

Technical Information Liquiphant FTL63 Density with Density Computer FML621

Vibronic



Density computer for liquids specifically for the food and life sciences industries

Application

The density measuring line can be used in liquid media and in hazardous areas. It is used for the following purposes:

- Density measurement
- Intelligent medium detection
- Reference density calculation
- To calculate the concentration of a liquid
- To convert values to different units such as °Brix, °Baumé, °API etc.

Advantages

- Certified, hygienic design (3-A, EHEDG)
- Proven conformity with materials standards, e.g. EC1935/2004, FDA, GB 4806, cGMP
- Measurement used directly in tanks or pipes without the need for additional pipework
- Integration of existing temperature measurements for temperature compensation
- Additional calculations, such as the concentration of a product, can be performed in the Density Computer FML621

Table of contents

About this document	4	Performance characteristics	28
Symbols	4	Reference operating conditions	28
Application	4	Measurement accuracy	28
Density measurement	4	Installation	29
Function and system design	7	Installation instructions for Liquiphant Density	29
Measuring principle	7	Density Computer FML621	32
System design	7	Environment	32
Specific density applications	7	Liquiphant Density	32
Measuring system	8	Density Computer FML621	33
Modularity	9	Liquiphant Density process	34
Electronic insert for density measurement	9	Process temperature range	34
Density Computer FML621	9	Thermal shock	34
Input of Liquiphant Density	9	Process pressure range	34
Measured variable	9	Pressure tightness	34
Measuring range	9	Solids contents	34
Output of Liquiphant Density	9	Mechanical construction	34
Output and input variants	9	Design, dimensions	34
Ex connection data	10	Dimensions	35
Input of Density Computer FML621	10	Weight	43
Measured variable	10	Materials	44
Measuring range	10	Surface roughness	45
Galvanic isolation	12	Mechanical construction of Density Computer	
Output of Density Computer FML621	12	FML621	45
Output signal	12	Terminal	45
Galvanic isolation	12	Dimensions	45
Current output, pulse output	12	Slots with extension cards	45
Switch output	13	Weight	46
Transmitter power supply and external power supply	13	Materials	46
Power supply of Liquiphant Density	13	User interface of Density Computer FML621	46
Terminal assignment	13	Display elements	46
Supply voltage	14	Operating elements	47
Power consumption	14	Remote operation	47
Current consumption	14	Real-time clock	47
Overvoltage protection	14	Certificates and approvals	47
Pulse signal in case of alarm	14	CE mark	47
Adjustment	15	Ex approval	47
Power supply of Density Computer FML621	15	Other standards and guidelines	47
Terminal assignment of density computer	15	Additional information on Liquiphant Density	47
Supply voltage	17	Ordering information	48
Power consumption	17	TAG	48
Power supply connection	17	Test reports, declarations and inspection certificates	49
Interface connection data	18	Accessories for Liquiphant Density	49
Slots, extension cards	19	Protective cover for aluminum dual compartment housing	49
Endress+Hauser-specific devices	19	Protective cover for aluminum single compartment housing	49
Connection of outputs	21	Weld-in adapter	50
Ethernet option	21	M12 socket	50
Extension cards (optional)	22		
Connecting the remote display and operating unit	27		

Accessories for Density Computer FML621 51
General 51
Extension cards 51
PROFINET® Interface 51

Documentation 52
Standard documentation 52
Supplementary device-dependent documentation 52

About this document

Symbols

Safety symbols



This symbol alerts you to a dangerous situation. Failure to avoid this situation will result in serious or fatal injury.



This symbol alerts you to a dangerous situation. Failure to avoid this situation can result in serious or fatal injury.




This symbol alerts you to a dangerous situation. Failure to avoid this situation can result in minor or medium injury.



This symbol contains information on procedures and other facts which do not result in personal injury.

Electrical symbols

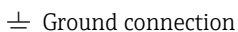


Grounded clamp, which is grounded via a grounding system.



Ground terminals, which must be grounded prior to establishing any other connections. The ground terminals are located on the inside and outside of the device.

Tool symbols

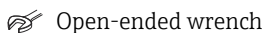

Symbols for certain types of information



Procedures, processes or actions that are permitted.



Procedures, processes or actions that are forbidden.



Indicates additional information



Reference to another section



1., 2., 3. Series of steps

Symbols in graphics

A, B, C ... View

1, 2, 3 ... Item numbers




Safe area (non-hazardous area)

Application

Density measurement

The Liquiphant Density measures the density of a liquid medium in pipes and tanks. The device is suitable for all Newtonian (purely viscous) fluids. In addition, the device is also suitable for use in hazardous areas.

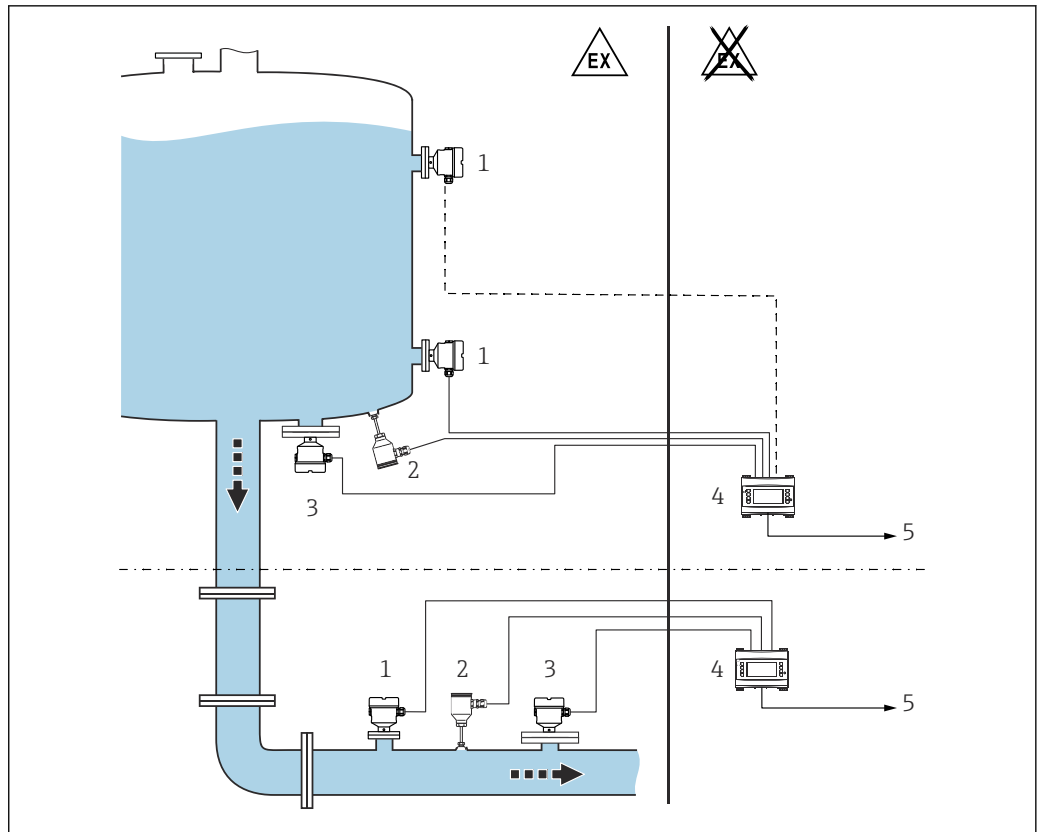

A0039632

1 Density measurement with FML621

- 1 Liquiphant Density → Pulse output
- 2 Temperature sensor, e.g. 4 to 20 mA output
- 3 Pressure transmitter 4 to 20 mA output required for changes in pressure >6 bar
- 4 Liquiphant Density Computer FML621 with display and operating unit
- 5 PLC



The measurement can be affected by:

- air bubbles at the sensor
- unit not fully covered by the medium
- solid media buildup on sensor
- high flow velocity in pipes
- severe turbulence in the pipe due to inlet and outlet runs that are too short
- corrosion of the fork
- non-Newtonian (not purely viscous) behavior of media

Application examples: basic unit

1 density measuring line, pressure and temperature-compensated

- 1 Liquiphant with FEL60D
- 1 temperature transmitter 4 to 20 mA
- 1 pressure transmitter 4 to 20 mA
- 1 output: density 4 to 20 mA
- 1 output: temperature 4 to 20 mA
- **Product structure:** FML621-xxxAAAxxxx
- **Number of inputs:** 4x pulse input, 0 to 20 mA or 4 to 20 mA
- **Number of outputs:** 1x SPST relay, 2x 0 to 20 mA or 4 to 20 mA

2 density measuring lines, temperature-compensated

- 2 Liquiphant with FEL60D
- 2 temperature transmitters 4 to 20 mA
- 1 output: density 4 to 20 mA
- 1 output: temperature 4 to 20 mA

- **Product structure:** FML621-xxxAAxxxx
- **Number of inputs:** 4x pulse input, 0 to 20 mA or 4 to 20 mA
- **Number of outputs:** 1x SPST relay, 2x 0 to 20 mA or 4 to 20 mA

Application examples: basic unit + 2 extension cards

3 density measuring lines, 2 x temperature-compensated, 1 x pressure- and temperature-compensated

- 3 Liquiphant with FEL60D
- 3 temperature transmitters 4 to 20 mA
- 1 pressure transmitter 4 to 20 mA
- 3 outputs: density 4 to 20 mA
- 3 outputs: temperature 4 to 20 mA
- 1 relay for medium detection
- **Product structure:** FML621-xxxBBxxxx
- **Number of inputs:** 8 x pulse input, 0 to 20 mA or 4 to 20 mA
- **Number of outputs:** 5 x SPST relay, 6x 0 to 20 mA or 4 to 20 mA

Application examples: medium detection

Distinguish between 2 media

- **Product structure:** FML621-xxxAAxxxx basic unit
- **Use of inputs:**
 - 1x FEL60D
 - 1x temperature 4 to 20 mA
- **Information content:**
 - 1 output: density 4 to 20 mA
 - 1 output: temperature 4 to 20 mA
 - 1 Relay



The medium detection can refer to concentrations or phase transitions

Distinguish between 3 media

- **Product structure:** FML621-xxxBAxxxx basic unit with additional relay card
- **Use of inputs:**
 - 1x FEL60D
 - 1x temperature 4 to 20 mA
- **Information content:**
 - 1 output: density 4 to 20 mA
 - 1 output: temperature 4 to 20 mA
 - 1 relay: display product 1
 - 1 relay: display product 2
 - 1 relay: display product 3



The relays can activate subsequent processes by triggering actuators.

Applications: density

Density measurement or concentration calculation with pump protection

- **Product structure:** FML621-xxxBAxxxx basic unit
- **Use of inputs:**
 - 1x FEL60D
 - 1x temperature 4 to 20 mA
- **Information content:**
 - 1 output: density 4 to 20 mA
 - 1 output: temperature 4 to 20 mA
 - 1 relay to switch off the pump




In addition to determining the density and concentration, pump protection can also be implemented by setting the appropriate switching frequency.

Function and system design

Measuring principle

A piezoelectric drive causes the tuning fork of the Liquiphant Density to vibrate at its resonance frequency. If the density of the liquid medium changes, the resonance frequency of the tuning fork also changes. The density of the medium has a direct impact on the resonance frequency of the tuning fork. As specific medium properties and mathematic relations are preprogrammed in the system, the Density Computer calculates the exact concentration of a medium.

System design

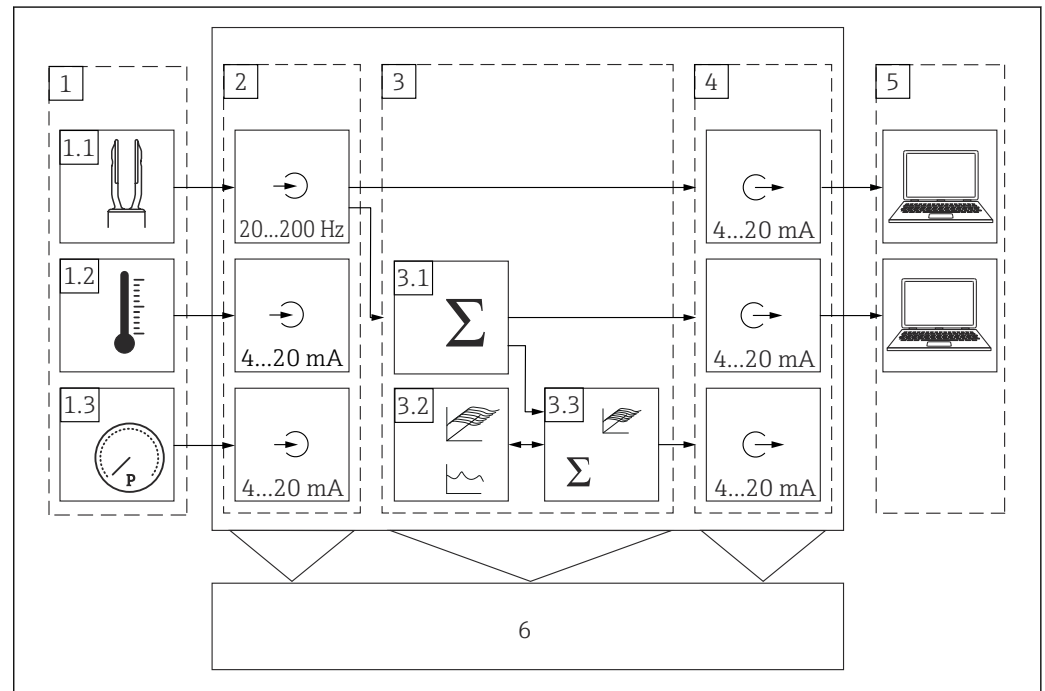

A0039647

2 Density Computer FML621, modular design diagram

- 1 External sensors
- 1.1 Liquiphant Density
- 1.2 Temperature sensor
- 1.3 Pressure sensor
- 2 Input modules, Density Computer FML621
- 3 Computing module, Density Computer FML621
- 3.1 Mathematical functions, e.g. density
- 3.2 2D, 3D curve
- 3.3 Mathematical functions, e.g. concentration, 3D linearization
- 4 Output modules, Density Computer FML621
- 5 Information processing - control room
- 6 Additional display

Specific density applications

The software modules available calculate the density from the frequency, temperature and pressure input variables.


Operating principle

The vibrating frequency of the tuning fork is reduced when the tuning fork is completely covered with liquid. Using additional information, such as temperature and pressure, it is possible to calculate the corresponding density of the medium. If the value by which the density has changed is known, the concentration of the medium can be determined using a function stored in the system. This value can be determined empirically or using existing tables, for example. The tables for converting density to concentration must be provided by the customer.

Additional software modules can calculate the density at reference temperature, compute the concentrations or detect media.

Reference density


In this module, the system refers to a reference temperature, such as 15 °C (59 °F) or 20 °C (68 °F). It must be known how the density of the medium changes at different temperatures.

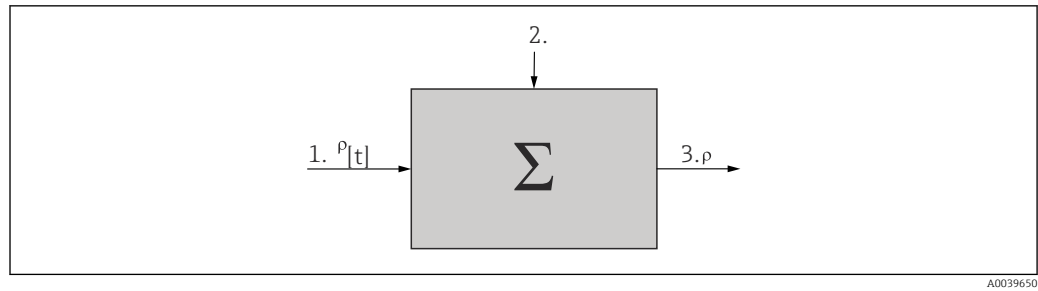

- 1 Input data: Table ρ [t]
- 2 Measured liquid medium: Temperature and density
- 3 Output: Calculated density ρ [standard]

Concentration


Using density and concentration curves already available or determined empirically, it is possible to determine the concentration when substances are continuously dissolved in a medium.

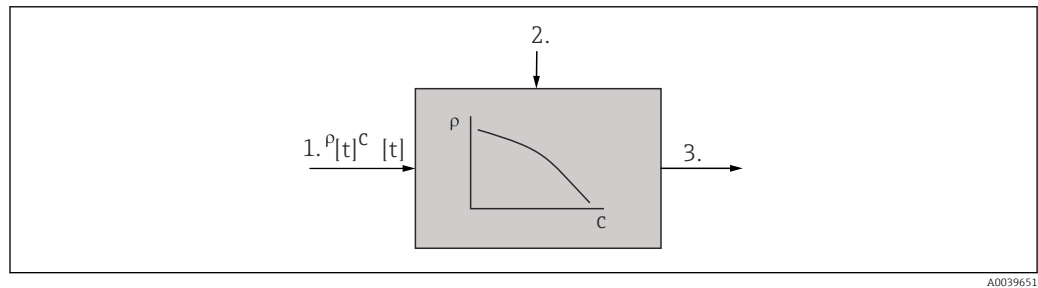

- 1 Input data: Table ρ, c [t]
- 2 Measured liquid medium: Temperature and density
- 3 Output: Calculated concentration

Medium detection

To be able to distinguish between two media, the density function - as a function of the temperature - can be stored for several media. This allows the system to differentiate between two media.

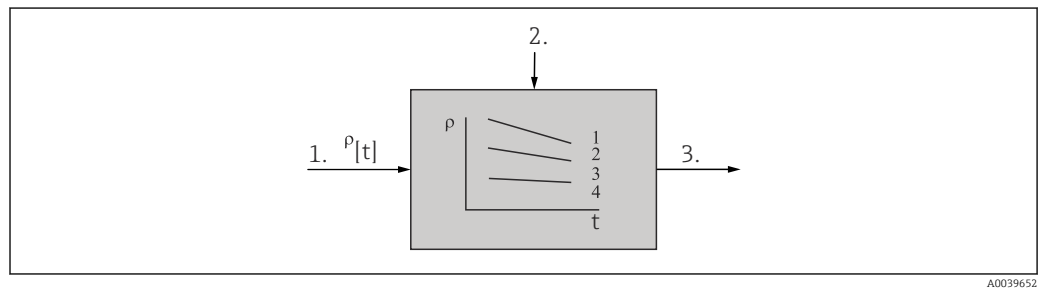

- 1 Input data: Tables ρ [t] for two liquid media
- 2 Measured liquid medium: Temperature and density
- 3 Output: Relay output, device

Measuring system

The Density Computer FML621 supplies power directly to connected two-wire transmitters. Intrinsically safe inputs and transmitter power supply units for current cards are optionally available for applications in hazardous areas. The inputs, outputs, limit values and display as well as commissioning and maintenance of the device are performed via a dot-matrix display with eight soft keys and a backlight as well as an RS232 or RS485 interface or ReadWin® 2000 PC software. In addition, it is possible to extend the device using additional extension cards.

A change in the background color indicates alarms or limit value violations. It is possible to configure the background color.

To use the telealarm function, we recommend common industrial modems that have an RS232 interface. The measured values and events or alarms are encoded and transmitted in accordance with the serial protocol. It is possible to query the protocol type.



The number of inputs, outputs, relays and transmitter power supply units contained in the basic unit can be individually extended using a maximum of three plug-in cards.

Modularity

- Measuring density of a liquid medium
- Liquiphant with electronic insert FEL60D and Density Computer FML621
- Also for hazardous areas
- Up to 5 density measuring lines can be operated with the Density Computer FML621. All slots must be fitted with plug-in cards.

Density Computer FML621 - specification**■ Input**

- FEL60D sensor
- 0 to 20 mA or 4 to 20 mA analog inputs
- 0 to 18 digital inputs
- 4 to 10 pulse inputs
- Temperature sensors (mA, mV, V, TC, RTD)

■ Output

- 2 to 8 Analog outputs 0 to 20 mA or 4 to 20 mA
- 2 to 8 Pulse outputs - active or passive
- 1 to 19 relays SPST, AC or DC

■ Communication

- Ethernet IP
- PSTN or GSM modem
- Serial bus RS232, RS485
- PROFIBUS® via coupler
- PROFINET® via coupler
- ReadWin® 2000 PC software

■ Power supply mode

- 4 - 10 devices, max. current consumption 30 mA
- 1 device, max. current consumption 80 mA

■ Internal memory

512 kB

■ Calculating functions

Pre-defined or editable

Electronic insert for density measurement

Electronic insert FEL60D

Density Computer FML621

Two-wire pulse output: Current pulses, superimposed on the power supply along the two-wire cabling

Input of Liquiphant Density

Measured variable

Density of liquids

Measuring range

Density range: 0.3 to 2 g/cm³ (18.7 to 125 lb/ft³) (0.3 to 2 SGU)

Output of Liquiphant Density

Output and input variants**2-wire density (FEL60D) for density measurement**

Connection to Density Computer FML621



For detailed information, see the Technical Information.

Ex connection data

See safety instructions (XA): All data relating to explosion protection are provided in separate Ex documentation and are available from the Downloads area of the Endress+Hauser website. The Ex documentation is supplied as standard with all devices approved for use in explosion hazardous areas.

Input of Density Computer FML621

Measured variable

- Voltage (analog and digital input)
- Current (analog input)
- PFM
- Pulse input

The following measured variables are implemented as an analog signal or pulse signal:

- Flow
- Level
- Pressure
- Temperature
- Density



Only Endress+Hauser flow sensors can be connected to the PFM input.

Not suitable for level and pressure measuring instruments.

Measuring range**Current input**

- 0 to 20 mA or 4 to 20 mA +10 % overrange
- Max. input current: 150 mA
- Input impedance: <10 Ω
- Accuracy 0.1 % of full scale value
- Temperature drift: 0.04 % / K (0.022 % / °F)
- Signal damping first-order low-pass filter, adjustable filter constants 0 to 99 s
- Resolution: 13 bit

Current input (U-I-TC card with intrinsically safe inputs)

- 0 to 20 mA or 4 to 20 mA +10 % overrange
- Max. input current: 80 mA
- Input impedance: ≈10 Ω
- Accuracy: 0.1 % of full scale value
- Temperature drift: 0.01 % / K 0.01 % / K (0.0056 % / °F)

PFM/pulse input

- Frequency range: 0.01 to 18 kHz
- Signal level - with approx. 1.3 kΩ series resistor at max. 24 V voltage level:
 - Low: 2 to 7 mA
 - High: 13 to 19 mA
- Measurement method: Period length or frequency measurement
- Accuracy: 0.01 % of reading
- Temperature drift: 0.01 % Over entire temperature range

Voltage input (digital input)

- Voltage level:
 - Low: -3 to 5 V
 - High: 12 to 30 V (as per IEC 61131-2)
- Input current typically: 3 mA with overload and reverse polarity protection
- Sampling frequency:
 - 4x4 Hz
 - 2x 20 kHz or 2x 4 Hz

Voltage input (analog input)

- Voltage: 0 to 10 V, 0 to 5 V, ±10 V, inaccuracy ±0.1 % of measuring range, input impedance >400 kΩ
- Voltage: 0 to 100 mV, 0 to 1 V, ±1 V, ±100 mV, inaccuracy ±0.1 % of measuring range, input impedance >1 MΩ
- Temperature drift: 0.01 % / K (0.0056 % / °F)

Resistance thermometer Pt100 as per ITS 90

- Measuring range: -200 to 800 °C (-328 to 1 472 °F)
- Accuracy: 4-wire connection 0.03 % of full scale value
- Type of connection: 3-wire or 4-wire system
- Measuring current: 500 µA
- Resolution: 16 bit
- Temperature drift: 0.01 % / K (0.0056 % / °F)

Resistance thermometer Pt500 as per ITS 90

- Measuring range: -200 to 250 °C (-328 to 482 °F)
- Accuracy: 4-wire connection 0.1 % of full scale value
- Type of connection: 3-wire or 4-wire system
- Measuring current: 500 µA
- Resolution: 16 bit
- Temperature drift: 0.01 % / K (0.0056 % / °F)

Resistance thermometer Pt1000 as per ITS 90

- Measuring range: -200 to 250 °C (-328 to 482 °F)
- Accuracy: 4-wire connection 0.08 % of full scale value
- Type of connection: 3-wire or 4-wire system
- Measuring current: 500 µA
- Resolution: 16 bit
- Temperature drift: 0.01 % / K (0.0056 % / °F)

Thermocouples (TC)

- J (Fe-CuNi), IEC 584
 - Measuring range: -210 to 999.9 °C (-346 to 1 832 °F)
 - Accuracy: ± (0.15 % of measuring range +0.5 K) from -100 °C
± (0.15 % of measuring range +0.9 °F) from -148 °F
- K (NiCr-Ni), IEC 584
 - Measuring range: -200 to 1 372 °C (-328 to 2 502 °F)
 - Accuracy: ± (0.15 % of measuring range +0.5 K) from -130 °C
± (0.15 % of measuring range +0.9 °F) from -202 °F
- T (Cu-CuNi), IEC 584
 - Measuring range: -270 to 400 °C (-454 to 752 °F)
 - Accuracy: ± (0.15 % of measuring range +0.5 K) from -200 °C
± (0.15 % of measuring range +0.9 °F) from -382 °F
- N (NiCrSi-NiSi), IEC 584
 - Measuring range: -270 to 1 300 °C (-454 to 1 386 °F)
 - Accuracy: ± (0.15 % of measuring range +0.5 K) from -100 °C
± (0.15 % of measuring range +0.9 °F) from -148 °F
- B (Pt30Rh-Pt6Rh), IEC 584
 - Measuring range: 0 to 1 820 °C (32 to 3 308 °F)
 - Accuracy: ± (0.15 % of measuring range +1.5 K) from 600 °C
± (0.15 % of measuring range +2.7 °F) from 1 112 °F
- D (W3Re/W25Re), ASTM E 998
 - Measuring range: 0 to 2 315 °C (32 to 4 199 °F)
 - Accuracy: ± (0.15 % of measuring range +1.5 K) from 500 °C
± (0.15 % of measuring range +2.7 °F) from 932 °F
- C (W5Re/W26Re), ASTM E 998
 - Measuring range: 0 to 2 315 °C (32 to 4 199 °F)
 - Accuracy: ± (0.15 % of measuring range +1.5 K) from 500 °C
± (0.15 % of measuring range +2.7 °F) from 932 °F
- L (Fe-CuNi), DIN 43710, GOST
 - Measuring range: -200 to 900 °C (-328 to 1 652 °F)
 - Accuracy: ± (0.15 % of measuring range +0.5 K) from -100 °C
± (0.15 % of measuring range +0.9 °F) from -148 °F

- U (Cu-CuNi), DIN 43710
 - Measuring range: -200 to 600 °C (-328 to 1112 °F)
 - Accuracy: $\pm$ (0.15 % of measuring range +0.5 K) from -100 °C
 $\pm$ (0.15 % of measuring range +0.9 °F) from -148 °F
- S (Pt10Rh-Pt), IEC 584
 - Measuring range: 0 to 1768 °C (32 to 3214 °F)
 - Accuracy: $\pm$ (0.15 % of measuring range +3.5 K) for 0 to 100 °C
 $\pm$ (0.15 % of measuring range +1.5 K) from 100 to 1768 °C
 $\pm$ (0.15 % of measuring range +6.3 °F) for 0 to 212 °F
 $\pm$ (0.15 % of measuring range +2.7 °F) for 212 to 2314 °F
- R (Pt13Rh-Pt), IEC 584
 - Measuring range: -50 to 1768 °C (-58 to 3214 °F)
 - Accuracy: $\pm$ (0.15 % of measuring range +3.5 K) for 0 to 100 °C
 $\pm$ (0.15 % of measuring range +1.5 K) from 100 to 1768 °C
 $\pm$ (0.15 % of measuring range +6.3 °F) for 0 to 212 °F
 $\pm$ (0.15 % of measuring range +2.7 °F) for 212 to 2314 °F

Galvanic isolation

The inputs between the individual extension cards and the basic unit are galvanically isolated ($\rightarrow$ 12).



With digital inputs, all terminal blocks are galvanically isolated from one another.

Output of Density Computer FML621

Output signal

Current, pulse, transmitter power supply (MUS) and switching output

Galvanic isolation

- The signal inputs and outputs are galvanically isolated from the supply voltage.
Testing voltage: 2.3 kV
- All signal inputs and outputs are galvanically isolated from one another.
Testing voltage: 500 V



The specified insulation voltage is the AC testing voltage U_{eff} , which is applied between the connections. Basis for assessment: IEC 61010-1, protection class II, overvoltage category II.

Current output, pulse output

Current output

- 0 to 20 mA or 4 to 20 mA +10 % overrange, reversible
- Max. loop current: 22 mA- short-circuit current
- Max. load: 750 Ω at 20 mA
- Accuracy: 0.1 % of full scale value
- Temperature drift: 0.1 % /10 K (0.056 % / 10 °F) ambient temperature
- Output ripple: <10 mV at 500 Ω for frequencies <50 kHz
- Resolution: 13 bit
- Error signals: 3.6 mA or 21 mA limit value in accordance NAMUR NE 43 - adjustable

Pulse output

- Basic unit:
 - Frequency range: up to 12.5 kHz
 - Voltage level: 0 to 1 V low, 12 to 28 V high
 - Min. load: 1 k Ω
 - Pulse width: 0.04 to 1000 ms
- Extension cards - digital passive, open collector:
 - Frequency range: up to 12.5 kHz
 - $I_{\text{max}} = 200$ mA
 - $U_{\text{max}} = 24$ V $\pm$ 15 %
 - $U_{\text{low/max}} = 1.3$ V at 200 mA
 - Pulse width: 0.04 to 1000 ms

Number of outputs

- 2x 0 to 20 mA or 4 to 20 mA / pulse - in basic unit
- Ethernet option: No current output present in the basic unit
- Max. number:
 - 8x 0 to 20 mA or 4 to 20 mA / pulse - depends on the number of extension cards
 - 6 x digital passive - depends on the number of extension cards

Signal sources

All existing multifunctional inputs and results of mathematic calculations can be freely assigned to the outputs.

Switch output

Function

Limit relay switches in the operating modes: Minimum or maximum safety, gradient, alarm, frequency or pulse, device error.

Switching behavior

Binary, switches when the limit value is reached - potential-free NO contact.

Switching capacity

Max. 250 V_{AC} 3 A / 30 V_{DC} 3 A



Do not combine line voltage and protective extra low voltage for extension card relays.

Switching frequency

Maximum 5 Hz

Threshold

User-programmable

Hysteresis

0 to 99 %

Sig. Source

All available inputs and calculated variables can be allocated freely to the switching outputs.

Number of switching cycles

> 100,000

Scan rate

500 ms

Number

- 1 relay - in the basic unit
- Max. number: 19 relays - depends on the number and type of extension cards

Transmitter power supply and external power supply

Transmitter power supply, terminals 81/82 or 81/83 - optional power extension cards 181/182 or 181/183

- Max. output voltage: 24 V_{DC} ±15 %
- Impedance: <345 Ω
- Max. loop current: 22 mA (at U_{out} >16 V)

FML621 technical data:

- HART® communication is not impaired
- Number: 3 MUS in the basic unit
- Max. number: 10 - depends on the number and type of extension cards

Additional power supply terminals 91/92, e.g. external display:

- Supply voltage: 24 V_{DC} ±5 %
- Max. current: 80 mA, short-circuit proof
- Number: 1
- Source resistance: <10 Ω

Power supply of Liquiphant Density

NOTICE


Operation with other switching units is not permitted.

Destruction of electronic components.

- ▶ Do not install the FEL60D electronic insert in devices that were originally used as point level switches.

Terminal assignment

The output signal of the density sensor is based on pulse technology. With the aid of this signal, the fork frequency is constantly forwarded to the Density Computer FML621.

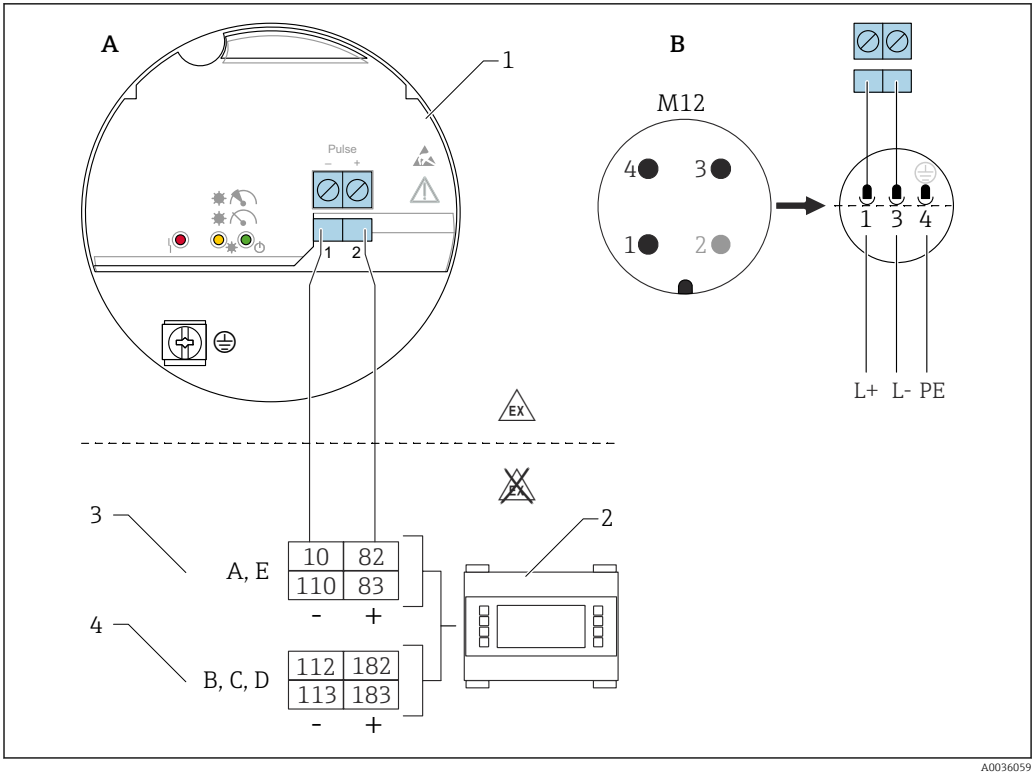

3 Connection diagram: electronic insert FEL60D connection to Density Computer FML621

- A Connection wiring with terminals
B Connection wiring with M12 plug in housing as per EN61131-2 standard
1 Electronic insert FEL60D
2 Density Computer FML621
3 Slots A, E with extension cards (already installed in the basic unit)
4 Slots B, C, D with extension cards (optional)

Supply voltage	U = 24 V _{DC} ±15 %, only suitable for connecting to the Density Computer FML621  The device must be powered by a voltage supply categorized as "CLASS 2" or "SELV".
Power consumption	P < 160 mW
Current consumption	I < 10 mA
Overvoltage protection	Overvoltage category I
Pulse signal in case of alarm	Output signal in case of power failure and damaged sensor: 0 Hz.

Adjustment

There are 3 different types of adjustment:

- **Standard adjustment (order configuration):**
Two fork parameters are determined at the factory to describe the sensor characteristics and are provided in the calibration report with the product. These parameters must be transmitted to the Density Computer FML621.
- **Special adjustment (select in the Product Configurator):**
Three fork parameters are determined at the factory to describe the sensor characteristics and are provided in the calibration report with the product. These parameters must be transmitted to the Density Computer FML621.
This type of adjustment achieves an even higher level of accuracy.
- **Field adjustment:**
With a field adjustment, the density determined by the user is transmitted to the FML621.



All the necessary parameters of the Liquiphant Density are documented in the **adjustment report** and in the **sensor pass**.


The documents are included in the scope of delivery.



Further information and the documentation currently available can be found on the Endress+Hauser website: www.endress.com → Downloads.

Power supply of Density Computer FML621


Terminal assignment of density computer

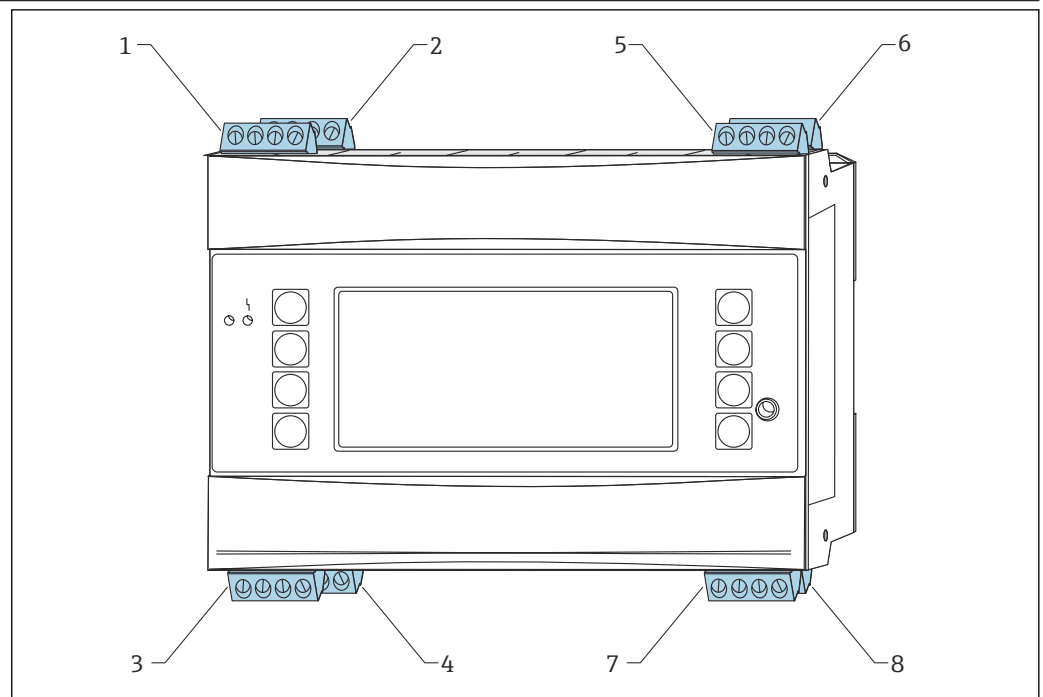

A0039654

4 Slot coding of basic unit

- 1 Slot A I - input
- 2 Slot A II - input
- 3 Slot A III - output
- 4 Slot A IV - output
- 5 Slot E I - input
- 6 Slot E II - input
- 7 Slot E III - output
- 8 Slot E IV - output

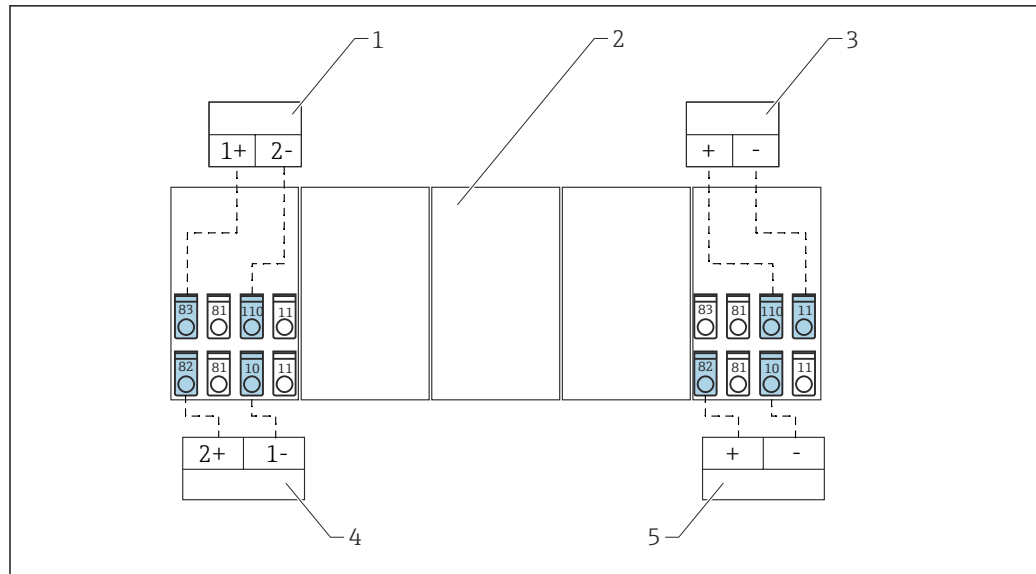


A0039655

5 Overview of connections - inputs

- 1 Passive sensor, e.g. pressure measurement
- 2 Slot for additional extension cards
- 3 Active sensor
- 4 Passive sensor, e.g. pressure measurement
- 5 Passive sensor, e.g. passive temperature transmitter

i Active sensor: The passing on of temperature information from a PLC is an example of why an active sensor may be connected.

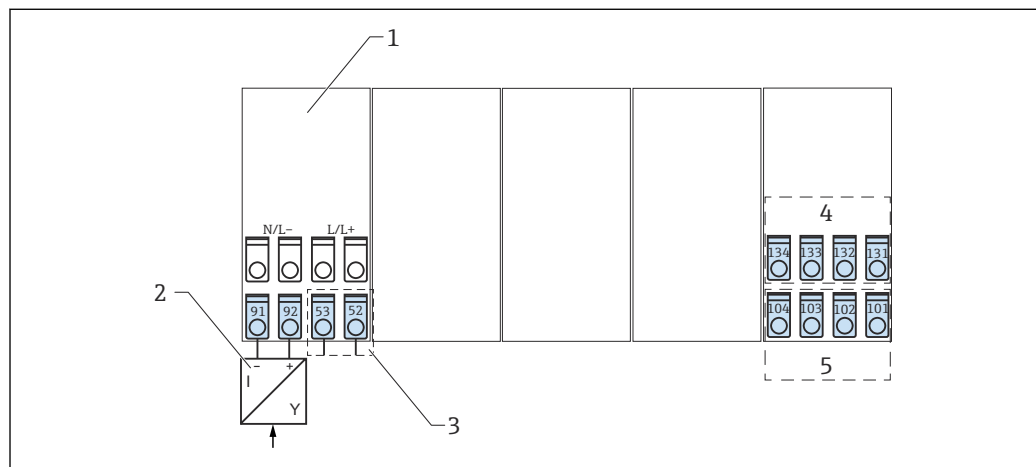

A0039656

6 Overview of connections - outputs

- 1 Extension card
- 2 Power supply for sensors
- 3 Relay contact
- 4 Pulse and current outputs - active
- 5 Bus interfaces

i With the Ethernet option, the current output or pulse output is not available at slot E.

Slot A I

Input: current or PFM or pulse input 1

- Terminal 10: (+) 0 to 20 mA or 4 to 20 mA, PFM, pulse input 1
- Terminal 11: ground for 0 to 20 mA or 4 to 20 mA, PFM, pulse input
- Terminal 81: ground, sensor power supply 1
- Terminal 82: 24 V sensor power supply 1

Slot A II

Input: current or PFM or pulse input 2

- Terminal 110: (+) 0 to 20 mA or 4 to 20 mA, PFM, pulse input 2
- Terminal 11: ground for 0 to 20 mA or 4 to 20 mA, PFM, pulse input
- Terminal 81: ground, sensor power supply 2
- Terminal 83: 24 V sensor power supply 2

Slot A III

Output: relay or additional sensor power supply

- Terminal 52: common relay (COM)
- Terminal 53: NO contact relay (NO)
- Terminal 91: ground, sensor power supply
- Terminal 93: +24 V sensor power supply

Slot A IV

Output: power supply

- Terminal L/L+: **L** for AC, **L+** for DC
- Terminal N/L-: **N** for AC, **L-** for DC

Slot E I

Input: current or PFM or pulse input 1

- Terminal 10: (+) 0 to 20 mA or 4 to 20 mA, PFM, pulse input 3
- Terminal 11: ground for 0 to 20 mA or 4 to 20 mA, PFM, pulse input
- Terminal 81: ground, sensor power supply 3
- Terminal 82: 24 V sensor power supply 3

Slot E II

Input: current or PFM or pulse input 2

- Terminal 110: (+) 0 to 20 mA or 4 to 20 mA, PFM, pulse input 4
- Terminal 11: ground for 0 to 20 mA or 4 to 20 mA, PFM, pulse input
- Terminal 81: ground, sensor power supply 4
- Terminal 83: 24 V sensor power supply 4

Slot E III

Output: RS485

- Terminal 101: (-) RxTx 1
- Terminal 102: (+) RxTx 1

Slot E III

Output: RS485 (optional)

- Terminal 103: (-) RxTx 2
- Terminal 104: (+) RxTx 2

Slot E IV

Output: current/pulse output 1

- Terminal 131: (-) 0/4 to 20 mA/pulse output 1
- Terminal 132: (+) 0/4 to 20 mA/pulse output 1

Slot E IV

 Ethernet, if the Ethernet option has been ordered.

Output: current/pulse output 2

- Terminal 133: (-) 0/4 to 20 mA/pulse output 2
- Terminal 134: (+) 0/4 to 20 mA/pulse output 2

 The inputs in the same slot are not galvanically isolated. There is a separation voltage of 500 V between the inputs and outputs in various slots. Terminals with identical second digits are bridged internally, for example terminals 11 and 81.

Supply voltage

- Low voltage power supply unit: 90 to 230 V_{AC} 50 to 60 Hz
- Extra-low voltage power supply unit: 20 to 36 V_{DC} or 20 to 28 V_{AC} 50 to 60 Hz

Power consumption

8 to 38 VA - depending on version and wiring.

Power supply connection

NOTICE

Destruction of electronic components.

- Check whether the supply voltage matches the voltage indicated on the nameplate of the device.

⚠ DANGER

Impermissible supply voltage

There is a high risk of physical injury and damage to the electronic components.


- For the device version with a supply voltage of 90 to 250 V, a switch must be installed in an easily accessible location. This switch is identified as a barrier in the power circuit of the device.

NOTICE

The power circuit of the device is not adequately protected.

Destruction of electronic components.

- Protect the power circuit with a 10-A fuse if the device is supplied with 90 to 250 V.

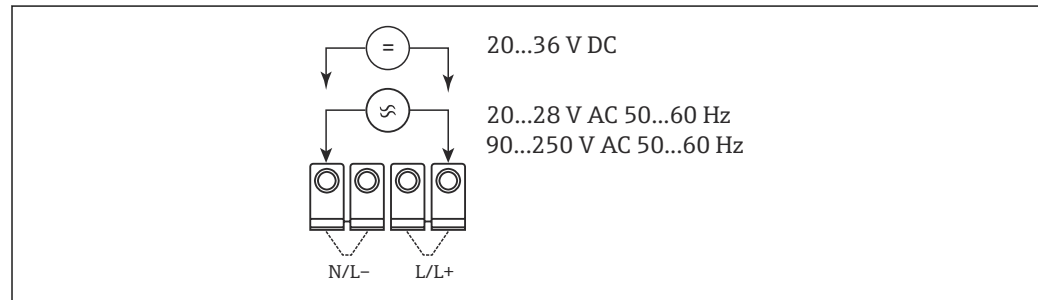

A0039657

 7 Power supply connection

Interface connection data

RS232

The RS232 interface is connected via an interface cable and a jack plug socket on the front of the housing.

- Connection: jack socket 3.5 mm (0.14 in), front
- Transmission protocol: ReadWin® 2000
- Transmission rate: max. 57 600 baud

RS485

- Connection: plug-in terminals 101 and 102
- Transmission protocol:
 - Serial: ReadWin® 2000
 - Parallel: open standard
- Transmission rate: max. 57 600 baud

PROFIBUS®, PROFINET®


- Optional connection of the Density Computer FML621 to PROFIBUS® or PROFINET® via the serial RS485 interface with the external HMS AnyBus protocol converter module for PROFIBUS® or PROFINET®
- Suitable protocol converter available as an accessory

Optional: additional RS485 interface

- Connection: plug-in terminals 103 and 104
- Transmission protocol and transmission rate as standard interface RS485

Optional: Ethernet interface

- Ethernet interface: 10/100 BaseT
- Connection type: RJ45
- Connection via shielded cable
- IP address output via Setup menu in device
- Connection to devices via an interface is only possible in office environments
- Safety distances: office device standard IEC 60950-1 must be taken into consideration
- Connection to a PC is possible via a "crossover" cable

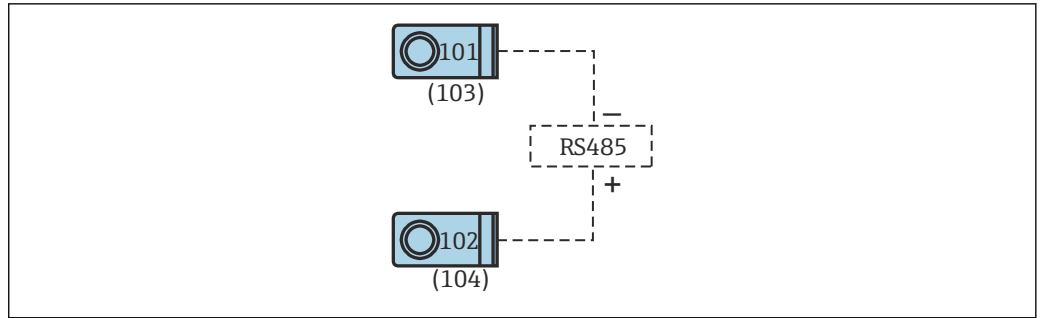

A0039688

8 Interface connection


Slots, extension cards

WARNING

The device is connected to the mains and is live.

Possibility of physical injury and destruction of electronic components.

- ▶ Ensure the device is in a de-energized state.
- ▶ Do not install or wire the device when it is connected to the mains voltage.

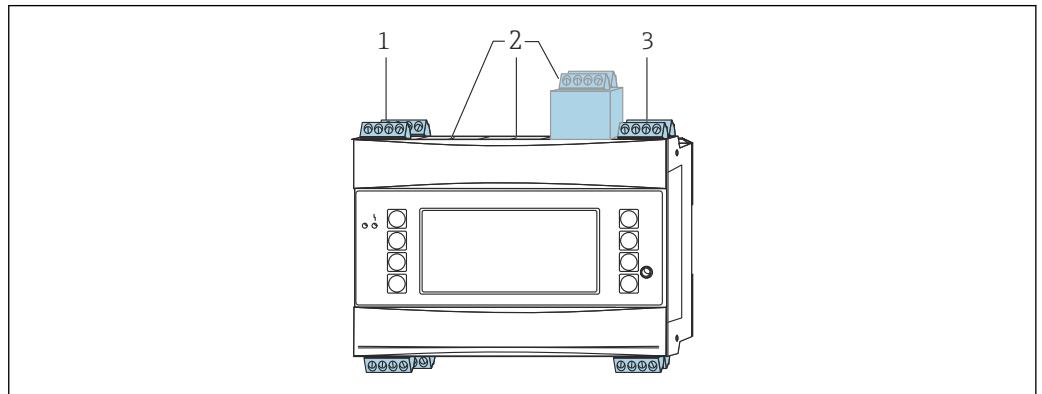

A0039653

9 Slots and extension cards in the density computer

- 1 Slot A, extension card already installed
- 2 Slots B, C, D can be extended with extension cards
- 3 Slot E, extension card already installed



The extension cards installed in slots A and E are an integral part of the basic unit.

Slots B, C and D can be expanded with additional extension cards.

Slots - specification

- Slot A:
 - Input: 2x density sensors, 0 to 20 mA or 4 to 20 mA
 - Output: 2x 0 to 20 mA or 4 to 20 mA
- Slots B, C, D:
 - Input: max. 10 analog inputs or 18 digital inputs
 - Output: max. 8 analog outputs or 6 digital outputs or 19 SPST relays
- Slot E:
 - Input: 2x density sensors 0 to 20 mA or 4 to 20 mA
 - Output: SPST relay

Endress+Hauser-specific devices




In the basic version of the Density Computer FML621, slots A and E are already fitted with extension cards.

Slots B, C and D can be expanded with additional extension cards.



The maximum cable length is 1 000 m (3 280.8 ft). The cable must be shielded to comply with EMC requirements. The maximum permitted power supply per core is 25 Ω.

Density sensor with a pulse output

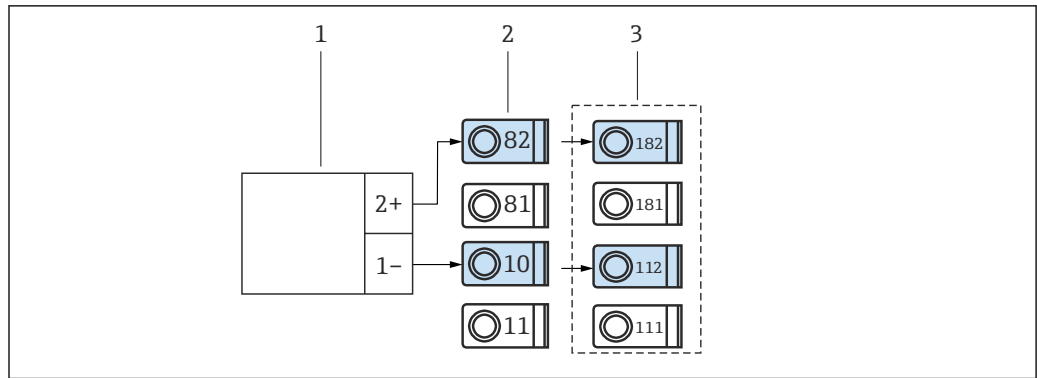


A0039671

10 Connecting the density sensor with pulse output

- 1 Density sensor
- 2 Slot A I
- 3 Additional slot B I

Temperature sensor via temperature head transmitter

i Connection of PT100, PT500 and PT1000 sensors is possible only via an optional extension card (in slot B, C, or D).

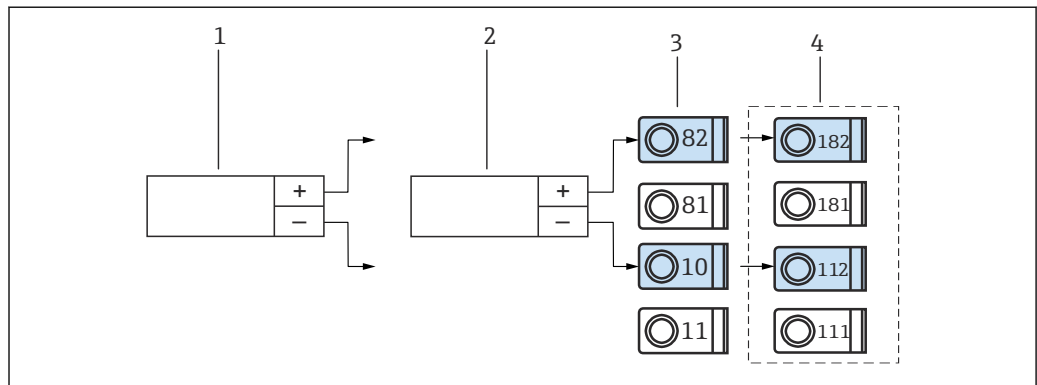


A0039673

11 Connecting the temperature sensor via the temperature head transmitter

- 1 Temperature transmitter 1
- 2 Temperature transmitter 2
- 3 Slot A I
- 4 Slot B I (optional extension card)

Pressure sensor with passive current output

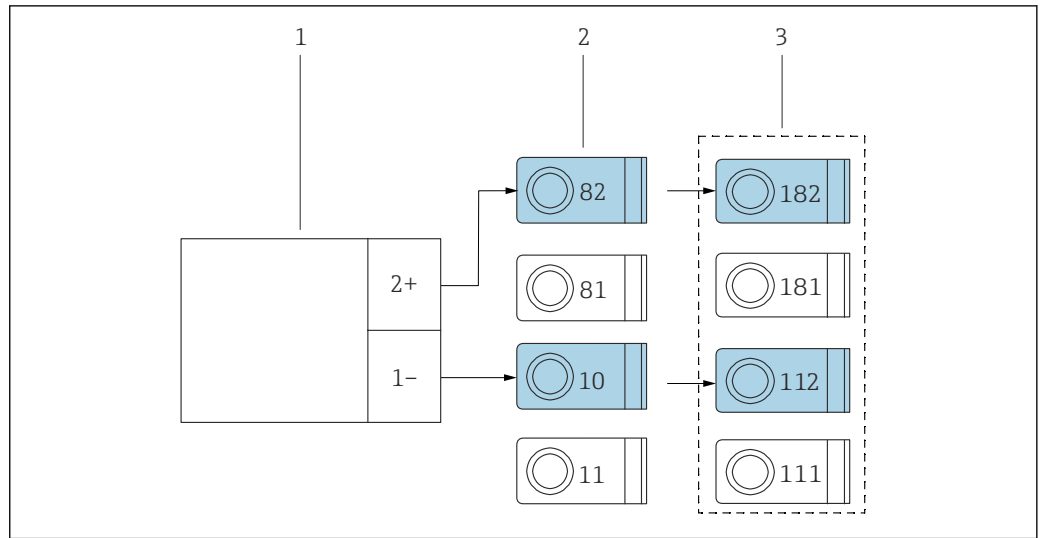


12 Connecting the pressure sensor with passive current output

- 1 Pressure transmitter
- 2 Slot A I
- 3 Slot B I (optional extension card)

Connection of outputs

The device has two galvanically isolated outputs or an Ethernet connection, which can be configured as an analog output or active pulse output. In addition, an output for connecting a relay and the option of transmitter power supply are available for each device. The number of outputs increases with the number of additionally installed extension cards (→ 22).

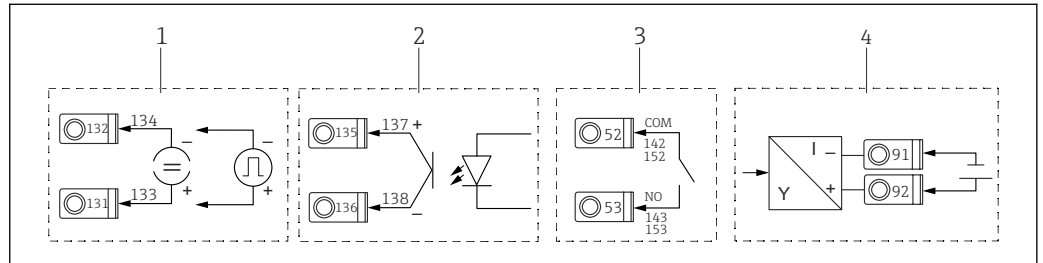

13 Connection of outputs


- 1 Active pulse and current outputs
- 2 Passive pulse output with open collector
- 3 Relay output (NO), e. g. slot A III
- 4 Transmitter power supply (MUS) output

Ethernet option

Ethernet connection

An IEEE 802.3-compatible connection on a shielded RJ45 connector on the underside of the device is available as a network connection. This can be used to connect the device to devices in the network environment with a hub or switch. For safety distances, the office equipment standard EN 60950 must be taken into account. The arrangement corresponds to a standard MDI interface (AT&T258), which means that a shielded 1:1 cable with a maximum length of 100 m (328 ft) can be used. The Ethernet interface is designed as a 10 and 100-BASE-T. Direct connection to a PC is possible with a crossover cable. Half-duplex and full-duplex data transmission are supported.

i If the Density Computer FML621 has an Ethernet interface, no analog outputs are available on the basic unit beyond slot E!

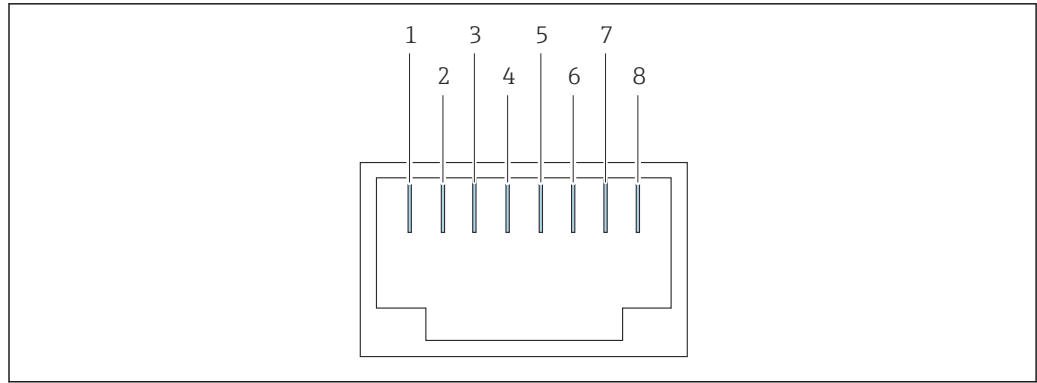

A0039690

14 RJ45 socket


- 1 Tx+
- 2 Tx-
- 3 Rx+
- 4 Not connected
- 5 Not connected
- 6 Rx-
- 7 Not connected
- 8 Not connected

LED indicators

Two LEDs below the plug-in connector indicate the status of the Ethernet interface:

- Yellow LED - connection signal
LED is lit when the device is connected to a network.
- Green LED - Tx/Rx
 - LED flashes when the device is sending or receiving data.
 - LED is continuously lit when the device is not sending or receiving data.

Extension cards (optional)

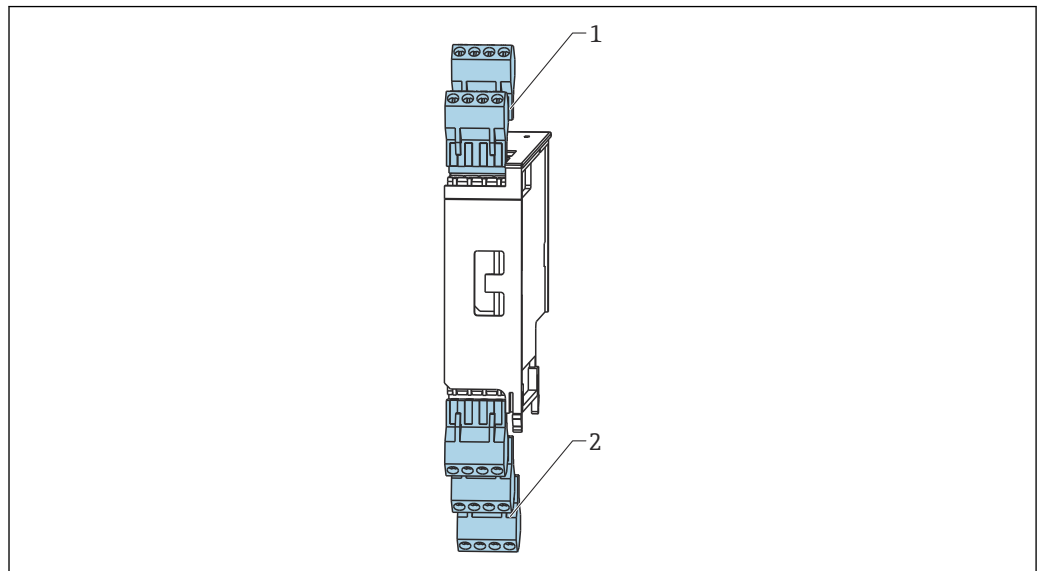

A0039691

15 Extension card with terminals (slots B, C and D)

- 1 Input: slots I, II
- 2 Output: slots III, IV, V

Terminal assignment, "Universal (FML621A-UA)" extension card with intrinsically safe inputs (FML621A-UB)

Slots B I, C I, D I

Input: current or PFM or pulse input 1

- Terminal 182: 24 V sensor power supply 1
- Terminal 112: (+) 0 to 20 mA or 4 to 20 mA, PFM, pulse input 1
- Terminal 111: ground for 0 to 20 mA or 4 to 20 mA, PFM, pulse input
- Terminal 181: sensor power supply ground 1

Slots B II, C II, D II

Input: current or PFM or pulse input 2

- Terminal 183: 24 V sensor power supply 2
- Terminal 181: sensor power supply ground 2
- Terminal 113: (+) 0 to 20 mA or 4 to 20 mA, PFM, pulse input 2
- Terminal 111: ground for 0 to 20 mA or 4 to 20 mA, PFM, pulse input

Slots B III, C III, D III

- Output: relay 1
 - Terminal 142: common relay (COM)
 - Terminal 143: NO contact relay (NO)
- Output: relay 2
 - Terminal 152: common relay (COM)
 - Terminal 153: NO contact relay (NO)

Slots B IV, C IV, D IV

Output: Current or pulse output - active

- Terminal 131: + 0 to 20 mA or 4 to 20 mA pulse output 1
- Terminal 132: - 0 to 20 mA or 4 to 20 mA pulse output 1
- Terminal 133: + 0 to 20 mA or 4 to 20 mA pulse output 2
- Terminal 134: - 0 to 20 mA or 4 to 20 mA pulse output 2

Slots B V, C V, D V

Output: Current or pulse output - passive


- Terminal 135: + pulse output 3 - open collector
- Terminal 136: - pulse output 3
- Terminal 137: + pulse output 4 - open collector
- Terminal 138: - pulse output 4

Terminal assignment, "Temperature (FML621A-TA)" extension card with intrinsically safe inputs (FML621A-TB)

Temperature sensors

Connection for Pt100, Pt500 and Pt1000.

 Terminals 116 and 117 must be bridged when connecting 3-wire sensors.

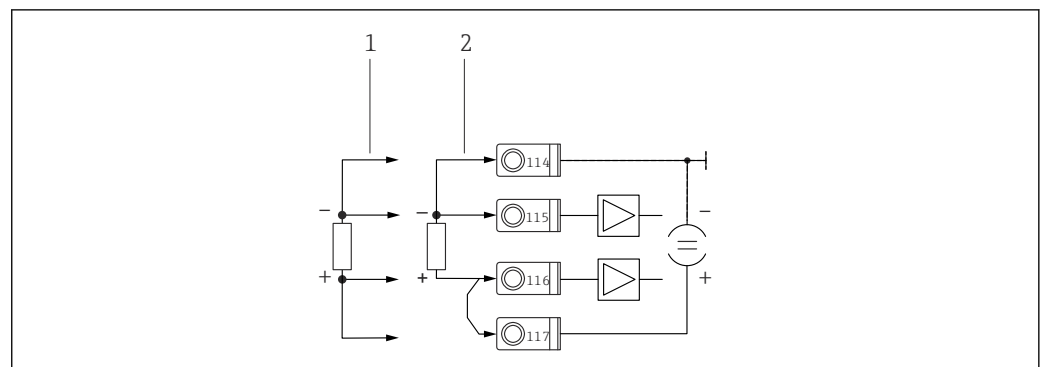

 16 Connecting the temperature sensor, optional temperature extension card e.g. in slot B (Slot B I)

- 1 4-wire input
- 2 3-wire input

Slots B I, C I, D I

Input: RTD input 1

- Terminal 117: + RTD power supply 1
- Terminal 116: + RTD sensor 1
- Terminal 115: - RTD sensor 1
- Terminal 114: - RTD power supply 1

Slots B II, C II, D II

Input: RTD input 2

- Terminal 121: + RTD power supply 1
- Terminal 120: + RTD sensor 1
- Terminal 119: - RTD sensor 1
- Terminal 118: - RTD power supply 1

Slots B III, C III, D III

- Output: relay 1
 - Terminal 142: common relay 1 (COM)
 - Terminal 143: NO contact relay 1 (NO)
- Output: relay 2
 - Terminal 152: common relay 2 (COM)
 - Terminal 153: NO contact relay 21 (NO)

Slots B IV, C IV, D IV

- Output: current or pulse output 1 - active
 - Terminal 131: + 0 to 20 mA or 4 to 20 mA
 - Terminal 132: - 0 to 20 mA or 4 to 20 mA
- Output: current or pulse output 2 - active
 - Terminal 133: + 0 to 20 mA or 4 to 20 mA
 - Terminal 134: - 0 to 20 mA or 4 to 20 mA


Slots B V, C V, D V

- Output: passive pulse output
 - Terminal 135: + pulse output 3 - open collector
 - Terminal 136: - pulse output 3
- Output: passive pulse output
 - Terminal 137: + pulse output 4 - open collector
 - Terminal 138: - pulse output 4

Terminal assignment of "Digital card (FML621A-DA)" extension card with intrinsically safe inputs (FML621A-DB)



The digital card has six intrinsically safe inputs. Terminals E1 and E4 can be defined as pulse inputs.

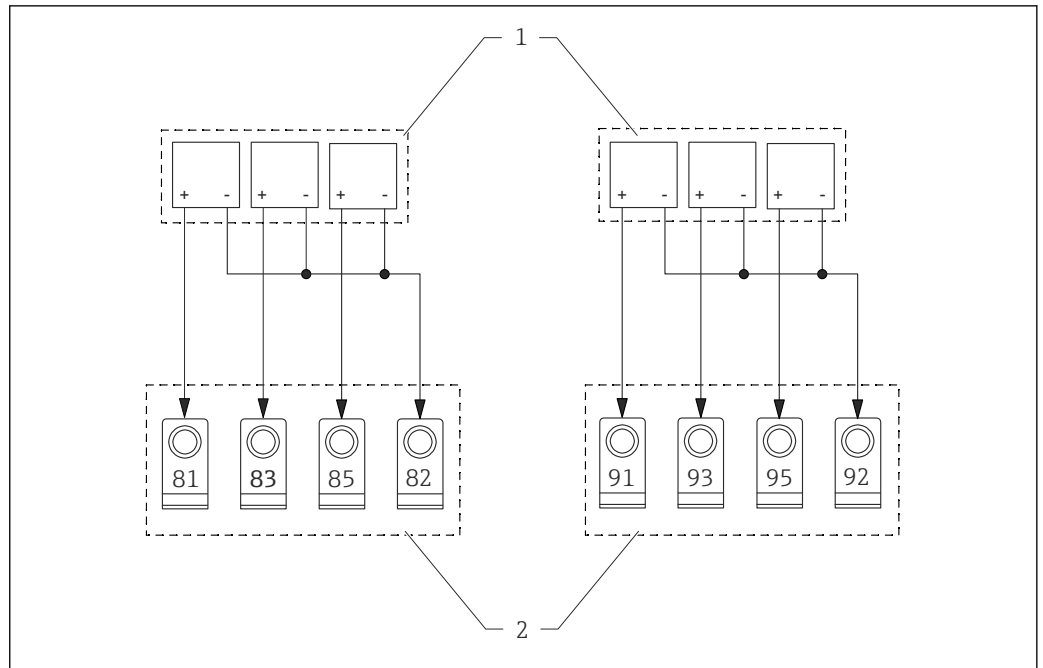

A0039694

17 Connecting the digital card

- 1 Digital input device
2 Terminal

i The current, PFM, pulse inputs or the RTD inputs in the same slot are not galvanically isolated. There is a separation voltage of 500 V between the aforementioned inputs and outputs in various slots.

Terminals with identical second digits are bridged internally.

Slots B I, C I, D I

Digital inputs E1 to 3

- Terminal 81: E1 20 kHz or 4 Hz as pulse input
- Terminal 83: E2 4 Hz
- Terminal 85: E3 4 Hz
- Terminal 82: signal ground E1 to 3

Slots B II, C II, D II

Digital inputs E4 to 6

- Terminal 91: E4 20 kHz or 4 Hz as pulse input
- Terminal 93: E5 4 Hz
- Terminal 95: E6 4 Hz
- Terminal 92: signal ground E4 to 6

Slots B III, C III, D III


- Output: relay 1
 - Terminal 142: common relay 1 (COM)
 - Terminal 143: NO contact relay 1 (NO)
- Output: relay 2
 - Terminal 152: common relay 2 (COM)
 - Terminal 153: NO contact relay 2 (NO)

Slots B IV, C IV, D IV

- Output: relay 3
 - Terminal 145: common relay 3 (COM)
 - Terminal 146: NO contact relay 3 (NO)
- Output: relay 4
 - Terminal 155: common relay 4 (COM)
 - Terminal 156: NO contact relay 4 (NO)

Slots B V, C V, D V

- Output: relay 5
 - Terminal 242: common relay 5 (COM)
 - Terminal 243: NO contact relay 5 (NO)
- Output: relay 6
 - Terminal 252: common relay 6 (COM)
 - Terminal 253: NO contact relay 6 (NO)

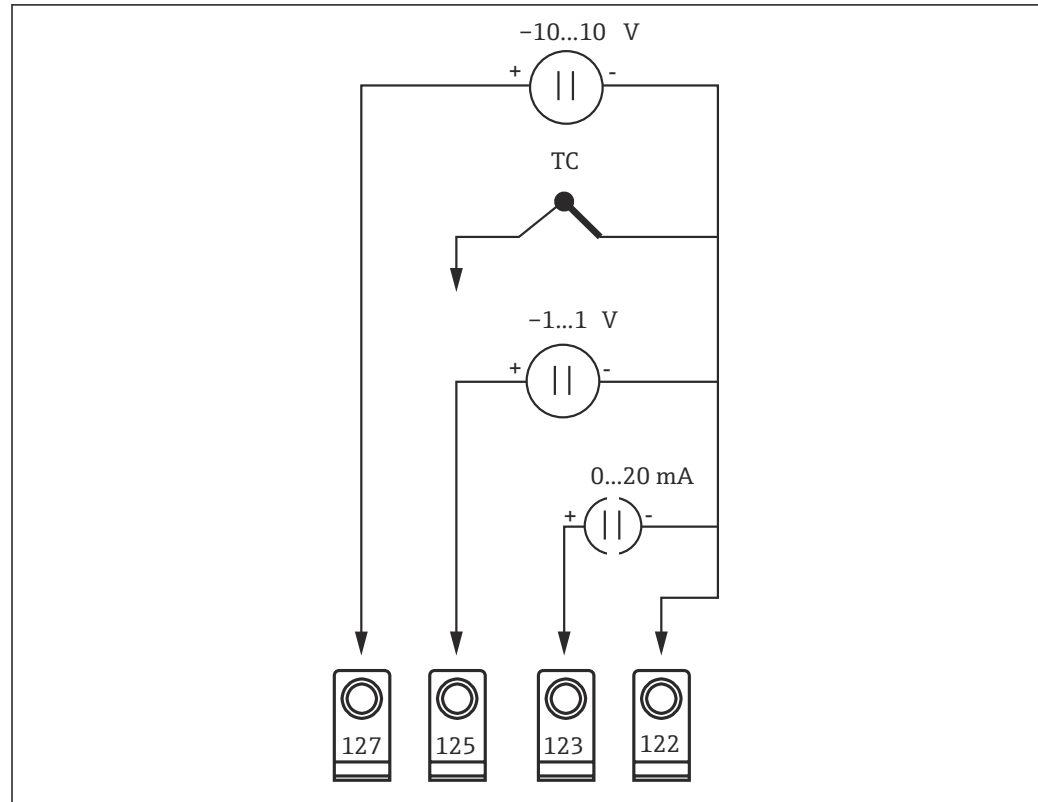Terminal assignment of extension card "U-I-TC card" with intrinsically safe inputs


A0039695

18 U-I-TC card



The card supports two input channels.

Channel 1 is supported by terminals 122, 123, 125 and 127.

Channel 2 is supported by terminals 222, 223, 225 and 227.

Slots B I, C I, D I
U-I-TC Input 1

- Terminal 127: -10 to +10 V input
- Terminal 125: -1 to +1 input, thermocouple
- Terminal 123: 0 to 20 mA input
- Terminal 122: input, signal ground

Slots B II, C II, D II
U-I-TC Input 2

- Terminal 227: -10 to +10 V input
- Terminal 225: -1 to +1 input, thermocouple
- Terminal 223: 0 to 20 mA input
- Terminal 222: input, signal ground

Slots B III, C III, D III

- Output: relay 1
 - Terminal 142: common relay 1 (COM)
 - Terminal 143: NO contact relay 1 (NO)
- Output: relay 2
 - Terminal 152: common relay 2 (COM)
 - Terminal 153: NO contact relay 2 (NO)

Slots B IV, C IV, D IV

- Output: current or pulse output 1 - active
 - Terminal 131: + 0 to 20 mA or 4 to 20 mA or pulse output 1
 - Terminal 132: - 0 to 20 mA or 4 to 20 mA pulse output 1
- Output: current or pulse output 2 - active
 - Terminal 133: + 0 to 20 mA or 4 to 20 mA or pulse output 2
 - Terminal 134: - 0 to 20 mA or 4 to 20 mA or pulse output 2

Slots B V, C V, D V

- Output: passive pulse output
 - Terminal 135: + pulse output 3 - open collector
 - Terminal 136: - pulse output 3
- Output: passive pulse output
 - Terminal 137: + pulse output 4 - open collector
 - Terminal 138: - pulse output 4

Connecting the remote display and operating unit

Description of functions

The remote display is an innovative addition to the powerful FML621 DIN rail device. The user has the opportunity to optimally install the arithmetic unit to suit the installation and mount the display and operating unit in a user-friendly way at easily accessible locations. The display can be connected to a DIN rail device both with and without an integrated display or operating unit. A 4-pin cable is supplied to connect the remote display with the basic unit. Other components are not necessary.



Please note the following:

- The remote display must be connected in order to use all of the functions of the operating unit
- Operating the unit solely with ReadWin® 2000 is not permitted
- Only ever connect one display or operating unit to the Density Computer FML621 (DIN rail device)

Installing the remote display or operating unit




The mounting location of the display unit must be free from vibrations.

The permitted ambient temperature during operation is -20 to +60 °C (-4 to +140 °F).

The device must be protected against high temperatures or heat.

Installing the display unit

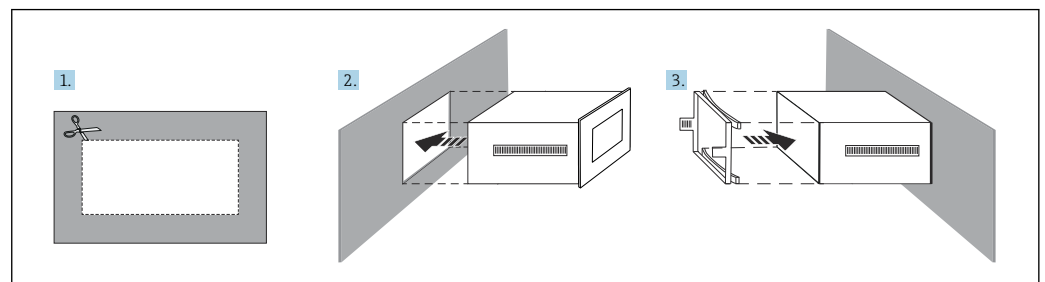


A0039697

19 Installing the display unit

1. Cut out a mounting opening with the following dimensions: 138 mm (5.43 in) x 68 mm (2.68 in), installation depth 43 mm (1.69 in).
2. Push the device, along with the sealing ring, through the cutout from the front.
3. Slide the securing frame over the rear of the housing and press it against the cabinet until the retaining clips click into place.
 - ↳ The display unit is now installed.

Wiring

The remote display and operating unit is connected directly to the basic unit using the enclosed cable.

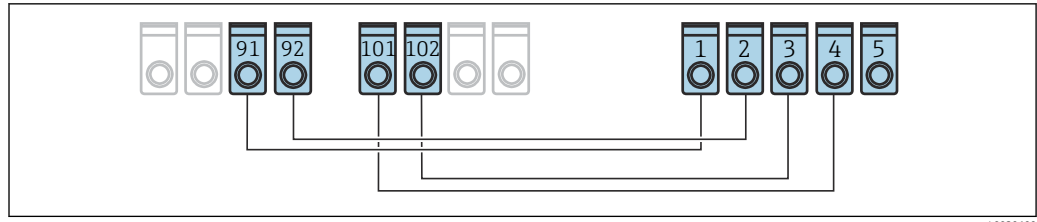

A0039699

20 Wire connections between remote display unit and basic unit.

- 1 Terminal GDN - remote display unit
- 2 Terminal 24 V_{DC} - remote display unit
- 3 Terminal + Rx Tx - remote display unit
- 4 Terminal - Rx Tx - remote display unit
- 5 Terminal PE - remote display unit
- 91 Terminal GND - slot A III - basic unit
- 92 Terminal 24 V_{DC} - slot A III - basic unit
- 101 Terminal - Rx Tx - slot E III - basic unit
- 102 Terminal + Rx Tx - slot E III - basic unit

Performance characteristics

Reference operating conditions

Normal operating conditions for special calibration and Liquiphant Density

- Medium: water H₂O
- Medium temperature: 0 to +80 °C (+32 to +176 °F), motionless fluid
- Ambient temperature: 24 °C (75 °F) ±5 °C (±9 °F)
- Moisture: max. 90 %
- Warm-up time: >30 min

Reference operating conditions of Density Computer FML621

- Power supply: 207 to 250 V_{AC} ±10 %, 50 Hz, ±0.5 Hz
- Warm-up time: >30 min
- Ambient temperature: +25 °C (+77 °F), ±5 °C (±9 °F)
- Humidity: 39 % ±10 % rF.

Measurement accuracy



The accuracy described here refers to the entire density measuring line.

General measuring conditions for accuracy data

- Measuring range: 0.3 to 2 g/cm³ (18.7 to 125 lb/ft³) (0.3 to 2 SGU)
- Observe the distance between the tuning fork and the surface of the medium (> 50 mm (1.97 in))
 see the "Orientation" section
- Measured error, temperature sensor: < 1 K
- Maximum viscosity: 350 mPa·s (3.5 P)
- Maximum flow velocity: 2 m/s (6.56 ft/s)
 - Laminar flow rate, bubble-free
 - For higher flow velocities, structural measures such as a bypass or an increase in pipe diameter must be taken to reduce the flow
- Process temperature: 0 to +80 °C (+32 to +176 °F) - validity of accuracy data
- Power supply as per FML621 specification
- Data as per DIN EN 61298-2
- Process pressure: -1 to +25 bar (-14.5 to +362.5 psi)

Measured error

1 g/cm³ (62.4 lb/ft³) = 1 SGU (Specific Gravity Unit)

- Standard adjustment: ±0.02 g/cm³ (±1.2 lb/ft³) (±1.2 % of the span 1.7 g/cm³ (106.1 lb/ft³), under general measuring conditions)
- Special adjustment: ±0.005 g/cm³ (±0.3 lb/ft³) (±0.3 % of span 1.7 g/cm³ (106.1 lb/ft³), under normal operating conditions)
- Field adjustment: ±0.002 g/cm³ (±0.1 lb/ft³), at operating point

Non-repeatability - reproducibility

1 g/cm³ (62.4 lb/ft³) = 1 SGU (Specific Gravity Unit)

- Standard adjustment: ±0.002 g/cm³ (±0.1 lb/ft³) (under general measuring conditions)
- Special adjustment: ±0.0007 g/cm³ (±0.04 lb/ft³) (under normal operating conditions)
- Field adjustment: ±0.002 g/cm³ (±0.1 lb/ft³), at operating point

Factors influencing accuracy data

-  Clean the sensor (CIP/SIP) if process temperatures of up to 140 °C (284 °F) persist over a long period
- All information relating to accuracy when determining the viscosity of liquids is based on Newtonian fluids
- A density measurement can be performed in the following liquids: gels, viscoelastic gels, non-Newtonian elastic fluids, pseudoelastic and plastic-viscous fluids.
- Long-term drift typically: ±0.00002 g/cm³ (±0.0012 lb/ft³) per day
- Temperature coefficient typically: ±0.0002 g/cm³ (±0.002 lb/ft³) per 10 K
- Flow velocity in pipes: > 2 m/s (6.56 ft/s)
- Buildup on fork
- Air bubbles in the case of vacuum applications or improper installation
- Incomplete coverage of the fork
- In the event of changes in pressure >6 bar (87 psi), a pressure measurement is necessary for compensation
- In the event of changes in temperature >1 K, a temperature measurement is necessary for compensation
- Mechanical stress, such as deformation of the tuning fork, can impair accuracy and must be avoided
- Devices exposed to mechanical stress must be replaced

Cyclic field calibration can take place depending on the accuracy required.

Installation


Installation instructions for Liquiphant Density



The following information is supplemented by additional documentation for the Liquiphant (Endress+Hauser website www.endress.com → Downloads)

Orientation

The mounting location must be selected such that the tuning fork and the membrane are always immersed in the medium.

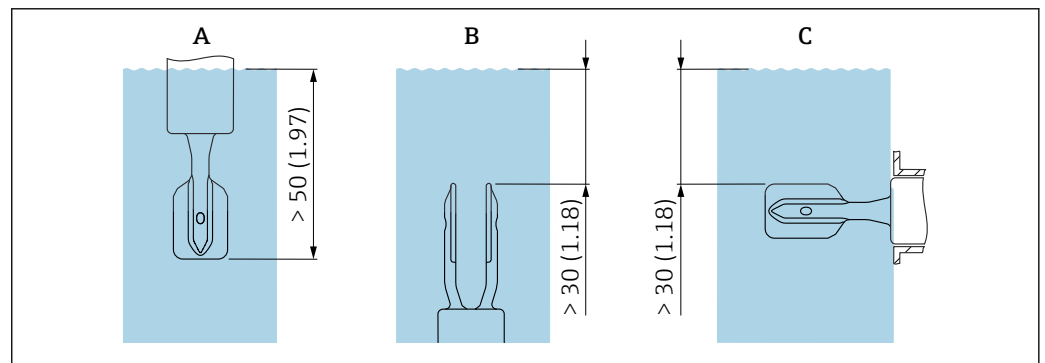

 21 Unit mm (in)

- A Installation from above
- B Installation from below
- C Installation from the side



- Avoid air bubbles in pipe or nozzle
- Ensure suitable ventilation

Input of correction factor "r"

The measurement result is affected if the distance between the tuning fork and the wall of the tank or pipe is very short:

- The medium has to flow around the tuning fork.
- The tuning fork of Liquiphant requires space to vibrate.

The measured error can be compensated for by entering a correction factor "r".

Pipe nominal diameters with internal measurements <44 mm (1.73 in) are not permitted!

 For detailed information, please refer to the relevant Operating Instructions.


Inlet and outlet runs

Inlet run

If possible, install the sensor as far upstream as possible, e.g. valves, T-pieces, elbows, flange elbows, etc.

To comply with the accuracy specification, the inlet run must meet the following requirements:

Inlet run: $\geq 5 \times \text{DN}$ (nominal diameter) - min. 750 mm (29.5 in)

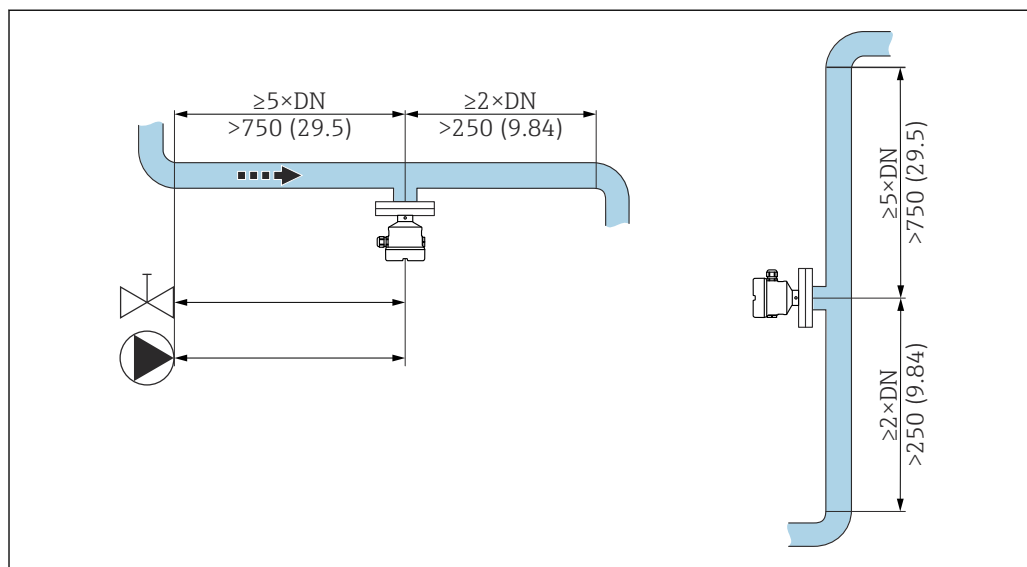


 22 Installing the inlet run. Unit of measurement mm (in)

Outlet run

To comply with the accuracy specification, the outlet run must meet the following requirements:

Outlet run: $\geq 2 \times \text{DN}$ (nominal diameter) - min. 250 mm (9.84 in)

The pressure and temperature sensor must be installed on the outlet side of the flow direction after the Liquiphant density sensor. When installing pressure and temperature measuring points downstream of the device, make sure the distance between the measuring point and the measuring device is sufficient.

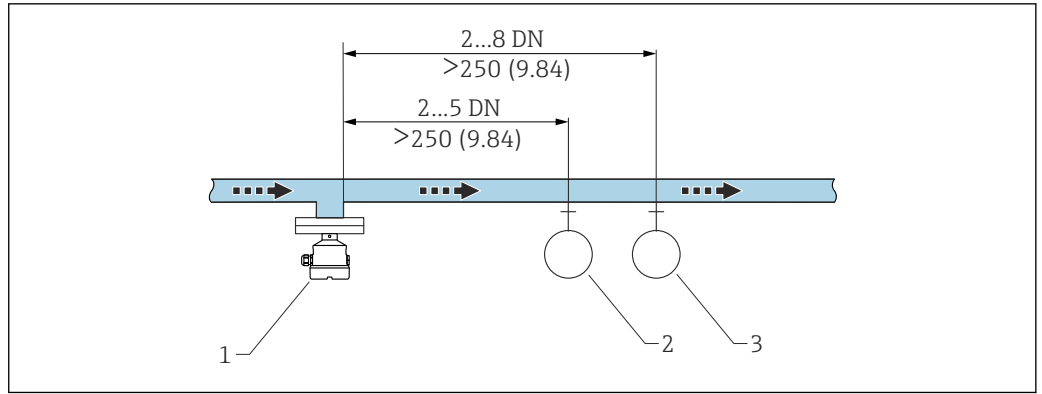

A0039701

23 Installing the outlet run. Unit of measurement mm (in)

- 1 Liquiphant density sensor
- 2 Pressure measuring point
- 3 Temperature measuring point


Installing the device in piping

NOTICE

Incorrect alignment of the tuning fork

Vortices and eddies can falsify the measuring result.

- ▶ Align the tuning fork in the flow direction for internal fixtures in pipes or tanks with an agitator.
- The flow velocity of the medium must not exceed 2 m/s (6.56 ft/s) during operation
- Flow velocity > 2 m/s: Separate the tuning fork from the direct flow of media by using structural features such as a bypass or pipe expansion to reduce the flow velocity to max. 2 m/s (6.56 ft/s)
- The flow will not be significantly impeded if the tuning fork is correctly aligned and the marking is pointing in the direction of flow.
- A marking on the process connection indicates the position of the tuning fork.
Threaded connection = dot on the hexagon head; flange = two lines on the flange.
The marking is visible when installed.

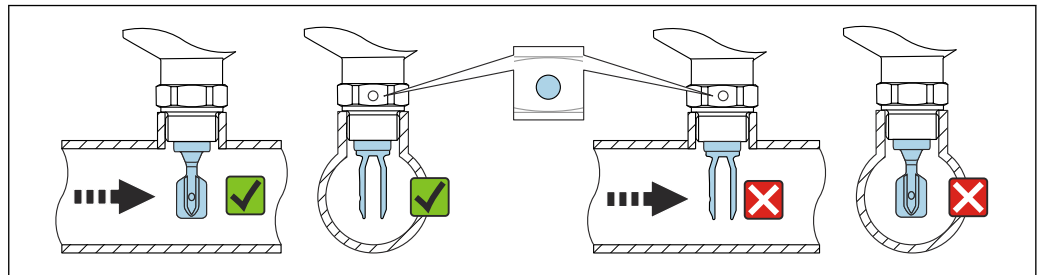

A0034851


24 Installation in pipes (take fork position and marking into consideration)

Aligning the cable entry

All housings can be aligned.

Housing without locking screw

The device housing can be rotated up to 350°.

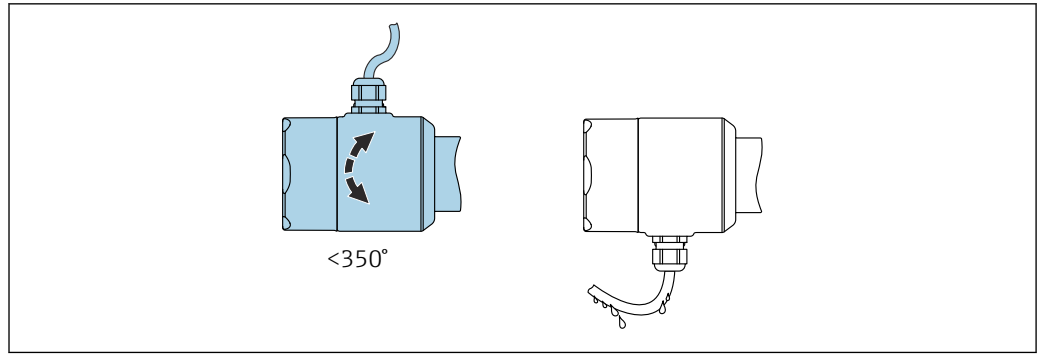

A0052359


25 Housing without locking screw with drip loop

Housing with locking screw



Housings with locking screw:

- The housing can be turned and the cable aligned by turning the locking screw.
- The locking screw is not tightened when the device is delivered.

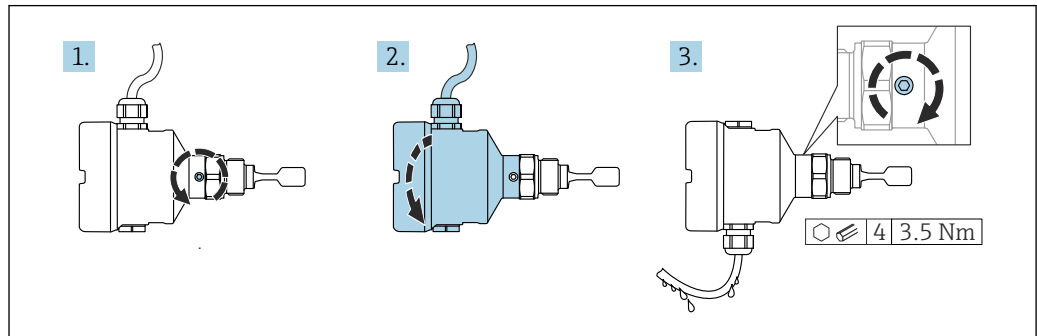

A0037347

26 Housing with external locking screw and drip loop

Density Computer FML621

Mounting location

Mount the device in a cabinet on a DIN rail as per IEC 60715.

Orientation

No restrictions.

Environment

Liquiphant Density

Ambient temperature range

–40 to 70 °C (–40 to 158 °F)

The minimum permitted ambient temperature of the plastic housing is limited to –20 °C (–4 °F); 'indoor use' applies for North America.

Outdoor operation in strong sunlight:

- Mount the device in a shaded location
- Avoid direct sunlight, particularly in warmer climatic regions
- Use a weather protection cover, can be ordered as an accessory



Further information on using the device in hazardous areas (ATEX) and on documentation that is currently available can be found on the Endress+Hauser website: www.endress.com → Downloads.

Humidity

Operation up to 100 %. Do not open in a condensing atmosphere.

Storage temperature

–40 to +80 °C (–40 to +176 °F)

Operating altitude

As per IEC 61010-1 Ed.3:

- Up to 2 000 m (6 600 ft) above sea level
- Can be extended to 3 000 m (9 800 ft) above sea level if overvoltage protection is used

Climate class

As per IEC 60068-2-38 test Z/AD

Degree of protection

Testing according to IEC 60529 and NEMA 250

IP68 test condition: 1.83 m H₂O for 24 h

Housing

See cable entries

Cable entries

- M20 coupling, plastic, IP66/68 NEMA Type 4X/6P
- M20 coupling, nickel-plated brass, IP66/68 NEMA Type 4X/6P
- M20 coupling, 316L, IP66/68 NEMA Type 4X/6P
- M20 coupling, 316L, hygienic, IP66/68/69 NEMA Type 4X/6P
- M20 thread, IP66/68 NEMA Type 4X/6P
- G ½ thread, NPT ½, NPT ¾ IP66/68 NEMA Type 4X/6P

Degree of protection for M12 plug

- When housing is closed and connecting cable is plugged in: IP66/67 NEMA Type 4X
- When housing is open or connecting cable is not plugged in: IP20, NEMA Type 1

NOTICE

M12 plug: Loss of IP protection class due to incorrect installation!

- The degree of protection only applies if the connecting cable used is plugged in and screwed tight.
- The degree of protection only applies if the connecting cable used is specified according to IP67 NEMA Type 4X.



If the "M12 plug" option is selected as the electrical connection, **IP66/67 NEMA Type 4X** applies for all housing types.

Pollution degree

Pollution degree 2

Density Computer FML621

Ambient temperature range

CAUTION

Extension cards generate additional heat.

Destruction of electronic components.

- Install additional ventilation with a minimum air current of 0.5 m/s (1.64 ft/s).

Temperature range: –20 to 50 °C (–4 to 122 °F).

Storage temperature

–30 to 70 °C (–22 to 158 °F)

Climate class

As per IEC 60654-1 Class B2/EN 1434 Class "C" - no condensation permitted.

Electrical safety

As per IEC 61010-1: operating environment at altitudes of < 2 000 m (6 560 ft) above sea level.

Degree of protection

- Basic unit: IP20
- Remote operating and display unit: front IP65

Electromagnetic compatibility**Interference emission**

IEC 61326 Class A

Interference immunity

- Power failure: 20 ms, no effect
- Starting current limitation: $I_{\max}/I_n < 50\%$ ($T 50\% \leq 50$ ms)
- Electromagnetic fields: 10 V/m (3.048 V/ft) as per IEC 61000-4-3
- Conducted HF: 0.15 to 80 Hz, 10 V as per IEC 61000-4-3
- Electrostatic discharge: 6 kV contact, indirect as per IEC 61000-4-2
 - Burst pulses - power supply: 2 kV as per IEC 61000-4-4
 - Burst pulses - signal: 1 kV/2 kV as per IEC 61000-4-4
 - Voltage peak - AC power supply: 1 kV/2 kV as per IEC 61000-4-5
 - Voltage peak - DC power supply: 1 kV/2 kV as per IEC 61000-4-5
 - Voltage peak - signal: 0.5 kV/1 kV as per IEC 61000-4-5

Liquiphant Density process

Process temperature range	0 to 80 °C (32 to 176 °F)
---------------------------	---------------------------

Thermal shock	≤ 120 K/s
---------------	----------------

Process pressure range	-1 to +25 bar (-14.5 to +362.5 psi)
------------------------	-------------------------------------

⚠ WARNING

The maximum pressure for the device depends on the lowest-rated element, with regard to pressure, of the selected component. This means that it is necessary to pay attention to the process connection as well as the sensor.

- ▶ Pressure specifications,  Technical Information, "Mechanical construction" section.
- ▶ Only operate the device within the specified limits!
- ▶ The Pressure Equipment Directive (2014/68/EU) uses the abbreviation "PS". The abbreviation "PS" corresponds to the MWP (maximum working pressure) of the device.

Pressure tightness	Up to vacuum
--------------------	--------------

Solids contents	$\varnothing \leq 5$ mm (0.2 in)
-----------------	----------------------------------

Mechanical construction


Design, dimensions**Device height**

The device height consists of the following components:

- Housing including cover
- Temperature spacer and/or pressure-tight feedthrough (second line of defense), optional
- Compact version, pipe extension or short pipe version
- Process connection

The individual heights of the components can be found in the following sections:

- Determine the height of the device and add the individual heights
- Take the installation clearance into consideration (space that is needed to install the device)

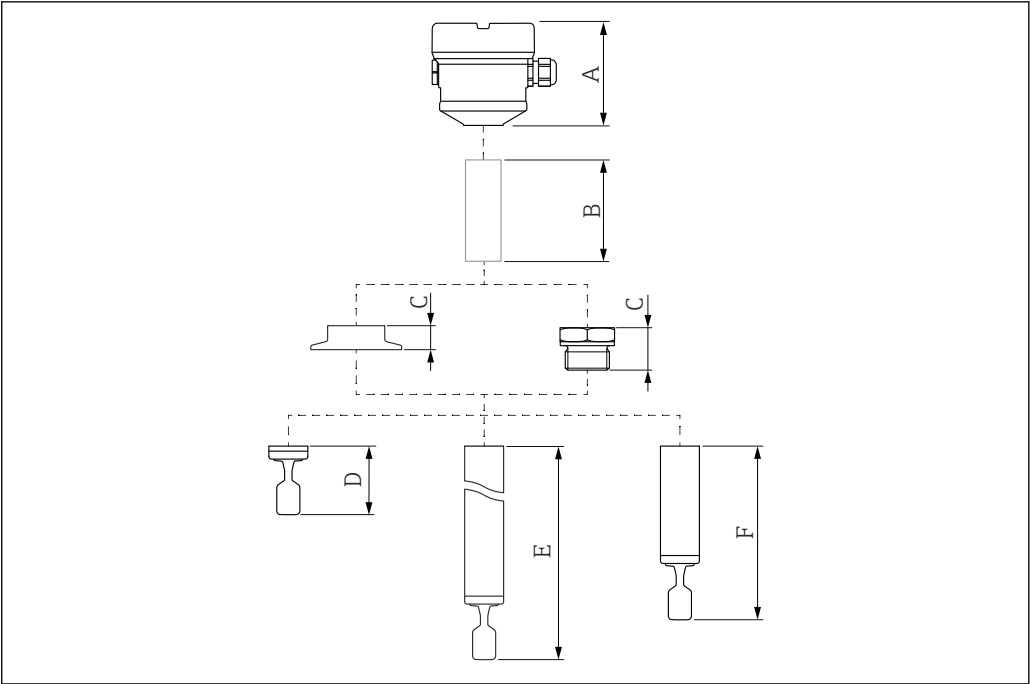

A0052410

27 Components to determine the device height


- A Housing including cover
- B Temperature spacer, pressure-tight feedthrough (optional)
- C Process connection
- D Probe design: compact version with tuning fork
- E Probe design: pipe extension with tuning fork
- F Probe design: short pipe version with tuning fork

Dimensions

Housing and cover

All housings can be aligned. The housing alignment can be fixed on housings with a locking screw.


Single compartment housing, plastic

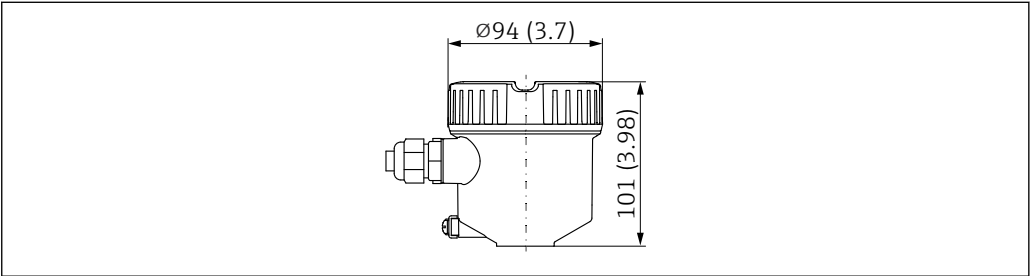

A0051909

28 Dimensions of single compartment housing, plastic; cover without sight glass. Unit of measurement mm (in)


Single compartment housing, aluminium, coated

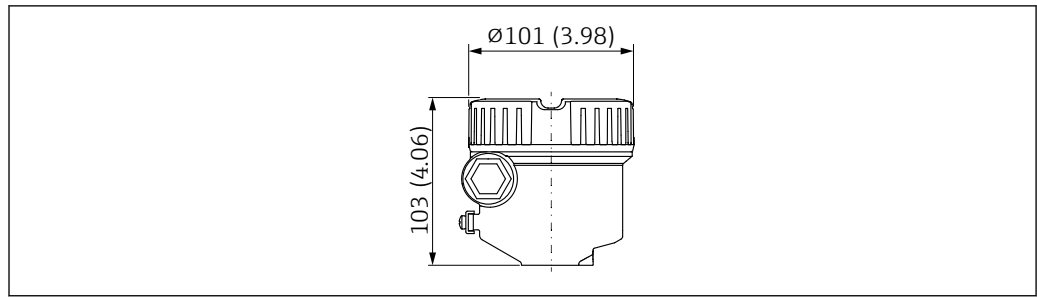

A0052195

- 29 Dimensions of single compartment housing, aluminium; cover without sight glass. Unit of measurement mm (in)

Single compartment housing, aluminum, coated (Ex d/XP, dust ignition-proof)

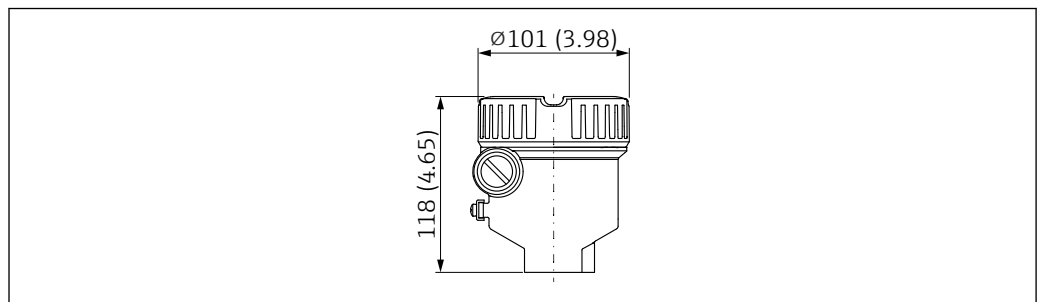


A0052194

- 30 Dimensions of single compartment housing, aluminum, coated; suitable for Ex d/XP, dust ignition-proof; cover without sight glass. Unit of measurement mm (in)

Single compartment housing, 316L, hygienic


- i** The housing with ground terminal and cover with cover lock are required for use in hazardous areas with a certain type of protection.

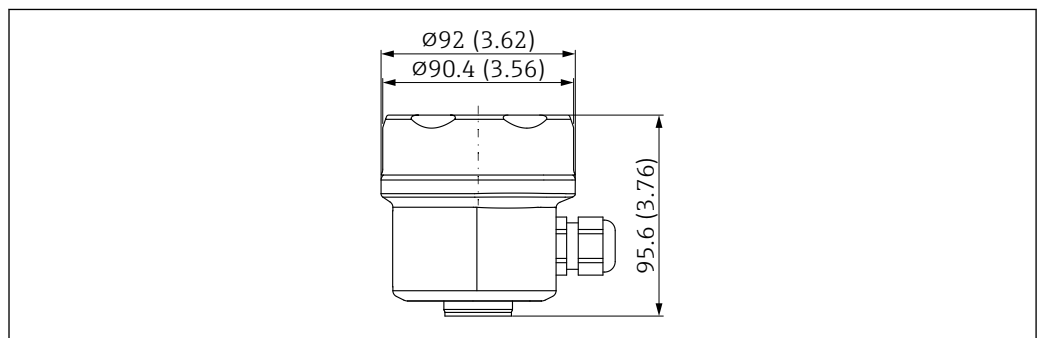

A0051667

- 31 Dimensions of single compartment housing, 316L, hygienic; cover without sight glass. Unit of measurement mm (in)

Dual compartment housing, L-shaped, aluminum, coated

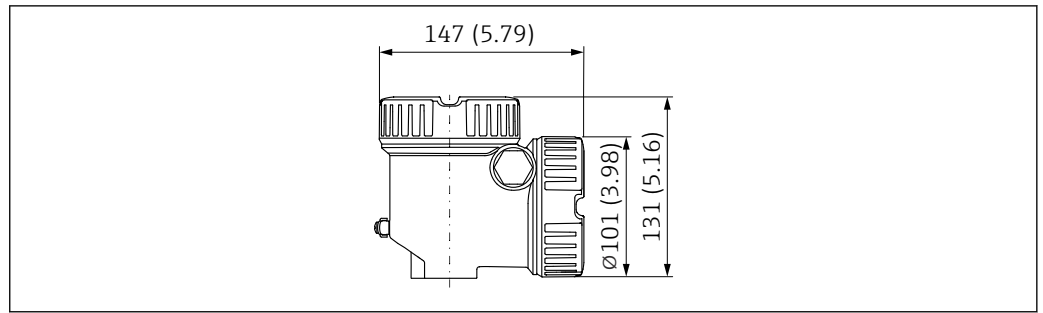

32 Dimensions of dual compartment housing, L-shaped, aluminum, coated; also with Ex d/XP, dust ignition-proof; cover without sight glass. Unit of measurement mm (in)

Ground terminal

- Ground terminal inside the housing, max. conductor cross-section 2.5 mm² (14 AWG)
- Ground terminal outside on the housing, max. conductor cross-section 4 mm² (12 AWG)

Cable glands

Cable outer diameter:

- Plastic: Ø5 to 10 mm (0.2 to 0.38 in)
- Nickel-plated brass: Ø7 to 10.5 mm (0.28 to 0.41 in)
- Stainless steel: Ø7 to 12 mm (0.28 to 0.47 in)
- Stainless steel, hygienic: Ø9 to 12 mm (0.35 to 0.47 in)




The scope of delivery comprises:

- 1 cable gland installed
- 1 cable gland sealed with dummy plug

Exceptions: For Ex d/XP, only threaded insertions are permitted.

Temperature spacer, pressure-tight feedthrough (optional)

The temperature spacer provides sealed insulation for the vessel.

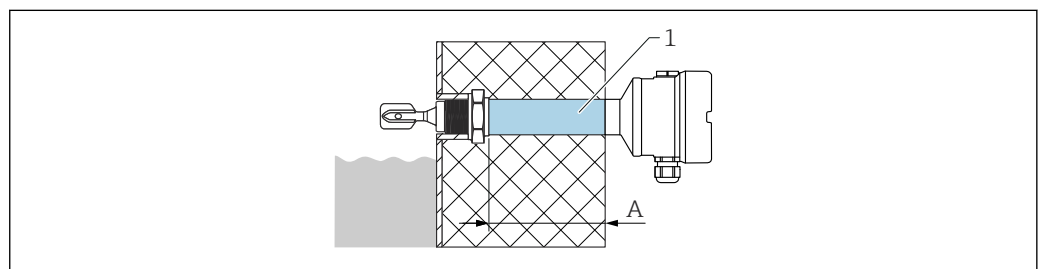

1 Temperature spacer and/or pressure-tight feedthrough with maximum insulation length
A 140 mm (5.51 in)

Product Configurator, feature "Sensor design":

- Temperature spacer
- Pressure-tight feedthrough (second line of defense)

If the sensor is damaged, this protects the housing from vessel pressures up to 100 bar (1 450 psi).



The "Pressure-tight feedthrough" version can only be selected in conjunction with the "Temperature spacer" option.


Probe design

Compact version

Sensor length L: depends on process connection



For more details, see the "Process connections" section.

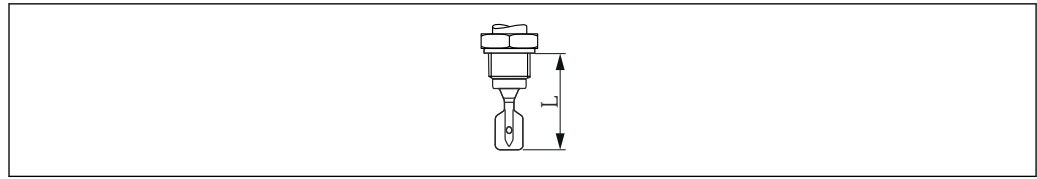

A0042435

33 Probe design: Compact version, sensor length L


Short pipe version

Sensor length L: depends on process connection

- Thread G 1 approx. 118 mm (4.65 in)
- Ingold, flush-mounted tank connection, DIN11851 DIN11864-1SMS1145 pipe union, DRD, Varivent, clamp/Tri-Clamp approx. 115 mm (4.53 in)
- Flush-mounted 1" (G 1 welding boss from Endress+Hauser): approx. 104 mm (4.09 in)

Pipe extension

- Sensor lengths L: 148 to 3 000 mm or 5.83 to 118.11 in
- Length tolerances L: < 1 m (3.3 ft) = -5 mm (-0.2 in), 1 to 3 m (3.3 to 9.8 ft) = -10 mm (-0.39 in)

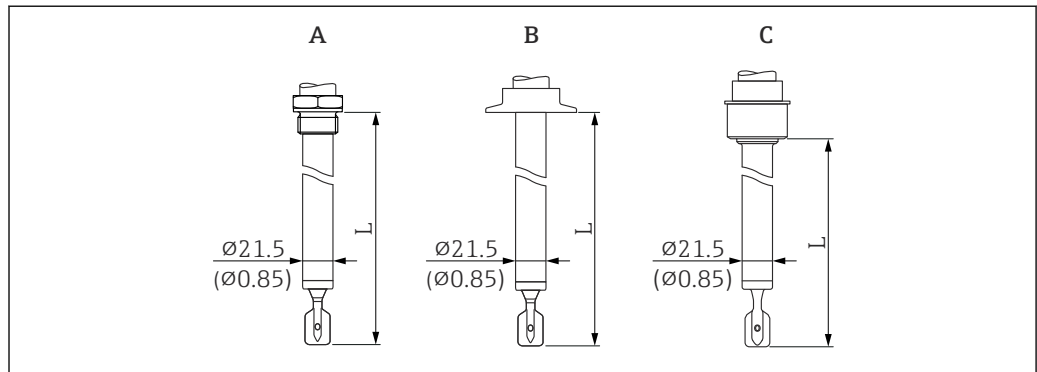

A0051989


34 Probe designs: pipe extension, short pipe version (sensor length L). Unit of measurement mm (in)

A Thread G 1

B E.g. clamp/Tri-Clamp, Varivent

C Flush-mounted tank connection for installation in weld-in adapter

Tuning fork

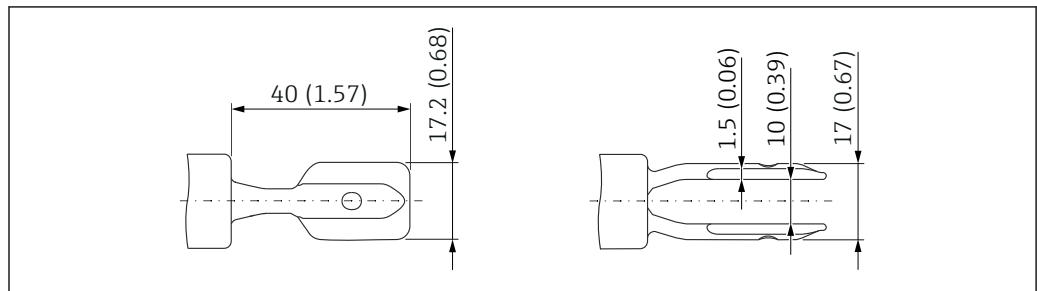

A0038269

35 Tuning fork. Unit of measurement mm (in)

Process connections


- i** Observe the permitted process temperature 0 to +80 °C (+32 to +176 °F). Higher temperatures affect the measurement accuracy.

Process connection, sealing surface

- Thread ISO228, G
- Ingold
- Flush-mounted tank connection

- DIN11851 pipe union
- DIN11864-1 pipe union
- DRD
- SMS1145 pipe union
- Varivent (Varinline)
- Clamp/Tri-Clamp

Height of process connection

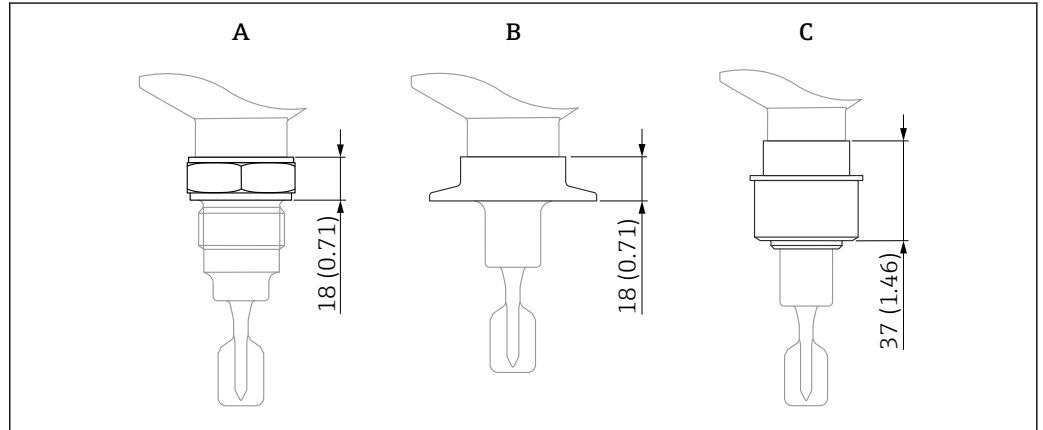

36 Maximum height specification for the process connections. Unit of measurement mm (in)


- A Process connection with threaded connection
 B For example: Clamp/Tri-Clamp, Varivent
 C Flush-mounted tank connection for installation in welding neck

Thread ISO228 G ¾ for installation in weld-in adapter

G ¾ with defined thread start for flush mounting in weld-in adapter

- Only for sensor design: compact version
- Material: 316L
- Pressure rating, temperature: ≤ 40 bar (580 psi), ≤ +100 °C (+212 °F)
- Pressure rating, temperature: ≤ 25 bar (363 psi), ≤ +150 °C (+302 °F)
- Weight: 0.2 kg (0.44 lb)
- Accessories: welding neck, optionally available as "Accessory enclosed"

i A seal is not included in the delivery. The maximum temperature and maximum pressure depend on the clamping ring and sealing element used (according to the process connection design). The lowest value applies in each case.

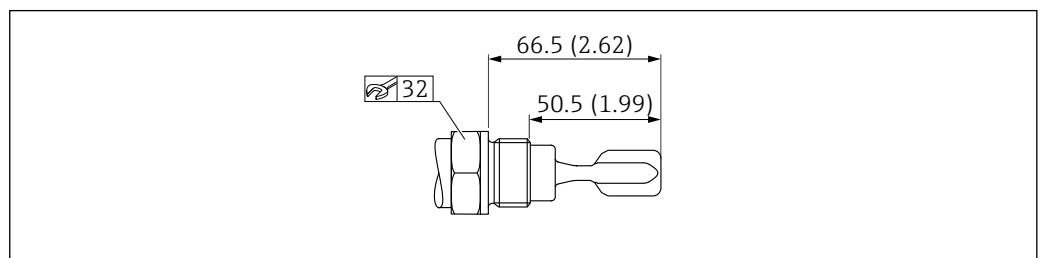


37 Thread ISO228 G ¾. Unit of measurement mm (in)

Thread ISO228 G 1 for installation in weld-in adapter

G 1 with defined thread start, comprising sealing surface for flush mounting in weld-in adapter

- Material: 316L
- Pressure rating, temperature: ≤ 40 bar (580 psi), $\leq +100$ °C (+212 °F)
- Pressure rating, temperature: ≤ 25 bar (363 psi), $\leq +150$ °C (+302 °F)
- Weight: 0.33 kg (0.73 lb)
- Accessories: welding neck, optionally available as "Accessory enclosed"

i A seal is not included in the delivery. The maximum temperature and maximum pressure depend on the clamping ring and sealing element used (according to the process connection design). The lowest value applies in each case.

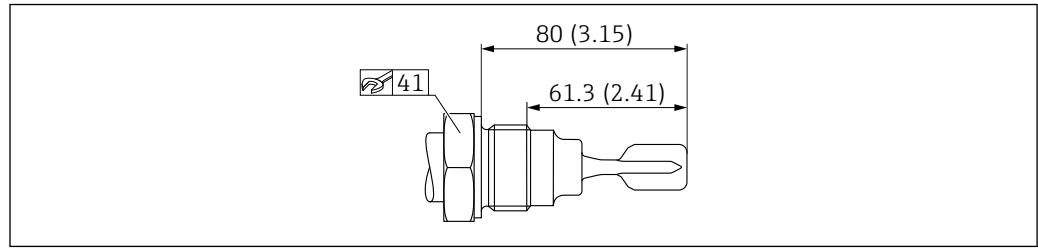


A0035551

38 Thread ISO228 G 1. Unit of measurement mm (in)

Ingold fitting

Ingold fitting 25 x 46 mm (2.52 in)

- Material: 316L
- Pressure rating: ≤ 16 bar (232 psi)
- Temperature: ≤ 150 °C (302 °F)
- Weight: 0.2 kg (0.44 lb)
- Scope of delivery: cap-nut G 1¼, seal

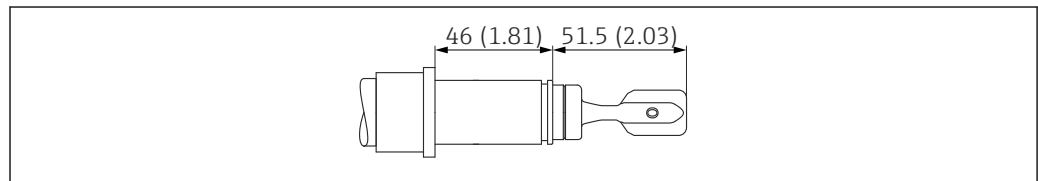


A0051991

39 Ingold fitting 25 x 46 mm (2.52 in). Unit of measurement mm (in)

Flush-mounted tank connection for installation in welding neck

- Material: 316L
- Pressure rating: ≤ 40 bar (580 psi) / ≤ 25 bar (363 psi)
- Temperature: ≤ 100 °C (212 °F) / ≤ 140 °C (284 °F)
- Weight: 0.44 kg (0.97 lb)
- Accessories: welding neck, optionally available as "Accessory enclosed"
- Scope of delivery: cap-nut, seal

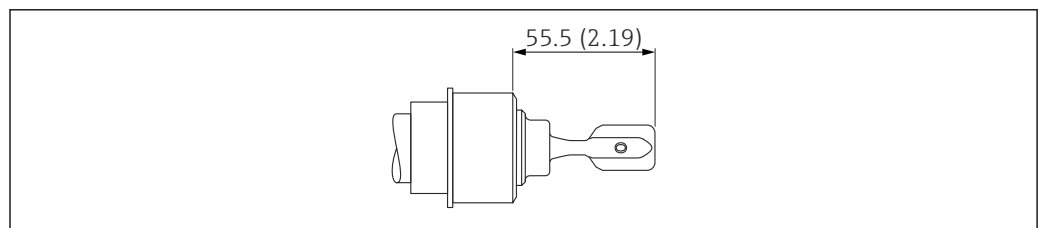

A0051993

40 Flush-mounted tank connection. Unit of measurement mm (in)

DIN11851 pipe union

DN32 PN25

- Material: 316L
- Slotted nut
- Pressure rating: ≤ 40 bar (580 psi) / ≤ 25 bar (363 psi)
- Temperature: ≤ 100 °C (212 °F) / ≤ 140 °C (284 °F)
- Weight: 0.3 kg (0.66 lb)


DN40 PN25

- Material: 316L
- Slotted nut
- Pressure rating: ≤ 40 bar (580 psi) / ≤ 25 bar (363 psi)
- Temperature: ≤ 100 °C (212 °F) / ≤ 140 °C (284 °F)
- Weight: 0.35 kg (0.77 lb)

DN50 PN25

- Material: 316L
- Slotted nut
- Pressure rating: ≤ 25 bar (363 psi)
- Temperature: ≤ 140 °C (284 °F)
- Weight: 0.47 kg (1.04 lb)

 A seal is not included in the delivery. The maximum temperature and maximum pressure depend on the clamping ring and sealing element used (according to the process connection design). The lowest value applies in each case.

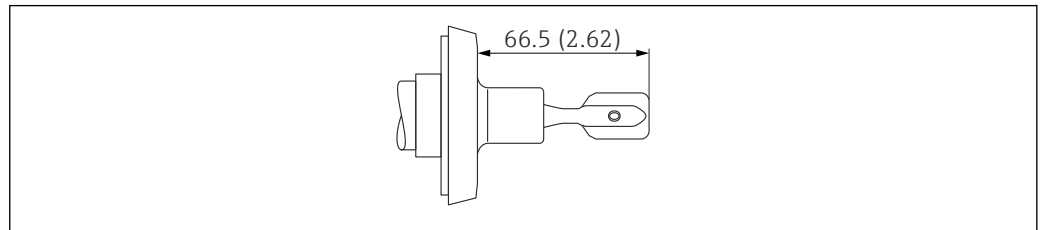

A0051995


 41 *DIN11851 pipe union. Unit of measurement mm (in)*

DIN11864-1 pipe union

DIN11864-1 A DN50 pipe DIN11850

- Material: 316L
- Slotted nut
- Pressure rating: ≤ 25 bar (363 psi)
- Temperature: ≤ 140 °C (284 °F)
- Weight: 0.47 kg (1.04 lb)

 A seal is not included in the delivery. The maximum temperature and maximum pressure depend on the clamping ring and sealing element used (according to the process connection design). The lowest value applies in each case.

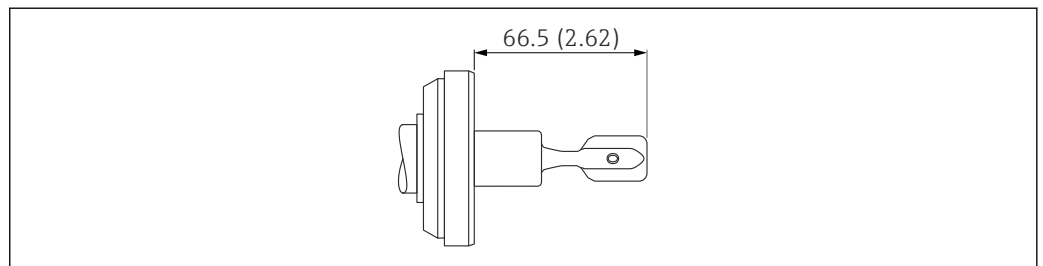

A0052381


 42 *DIN11864-1 pipe union. Unit of measurement mm (in)*

DRD

DRD 65 mm (2.56 in)

- Material: 316L
- Pressure rating: ≤ 40 bar (580 psi) / ≤ 25 bar (363 psi)
- Temperature: ≤ 100 °C (212 °F) / ≤ 140 °C (284 °F)
- Weight: 0.43 kg (0.95 lb)
- Accessories: welding flange with PTFE flat seal, optionally available as "Accessory enclosed"

i A seal is not included in the delivery. The maximum temperature and maximum pressure depend on the clamping ring and sealing element used (according to the process connection design). The lowest value applies in each case.

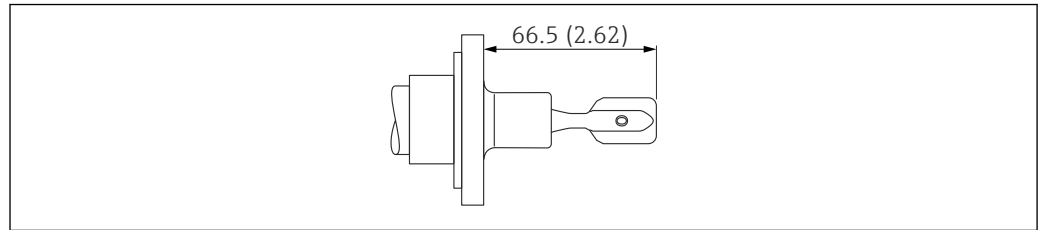

A0051992


43 DRD. Unit of measurement mm (in)

SMS1145 pipe union

SMS 2" PN25

- Material: 316L
- Pressure rating: ≤ 25 bar (363 psi)
- Temperature: ≤ 140 °C (284 °F)
- With cap-nut
- Weight: 0.33 kg (0.72 lb)

i A seal is not included in the delivery. The maximum temperature and maximum pressure depend on the clamping ring and sealing element used (according to the process connection design). The lowest value applies in each case.

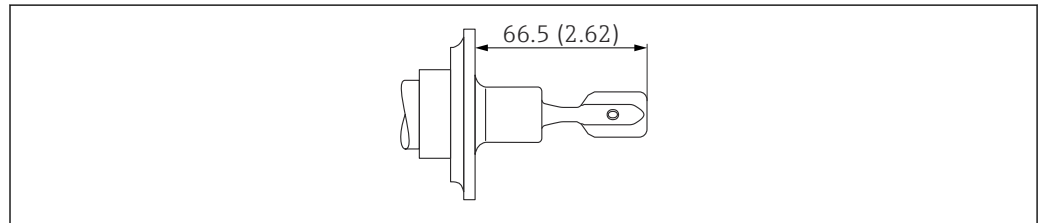

A0051994


44 SMS1145 pipe union. Unit of measurement mm (in)

Varivent (Varinline)

Varivent N pipe DN65-162 PN25

- Material: 316L
- Pressure rating: ≤ 25 bar (363 psi)
- Temperature: ≤ 150 °C (302 °F)
- Suitable for GEA Tuchenhausen
- Weight: 0.72 kg (1.59 lb)

i A seal is not included in the delivery. The maximum temperature and maximum pressure depend on the clamping ring and sealing element used (according to the process connection design). The lowest value applies in each case.

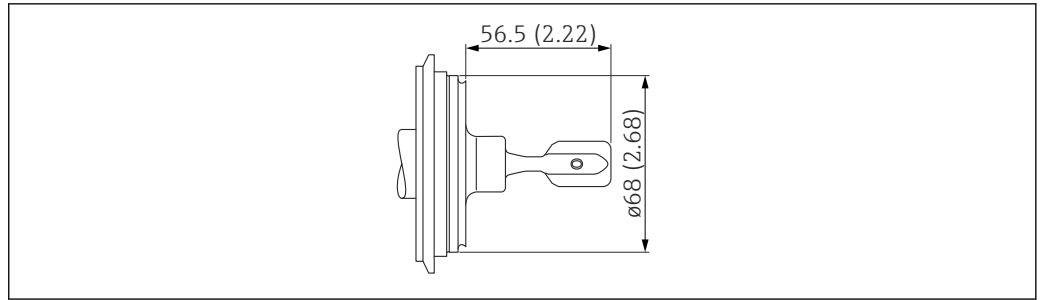

A0051996

45 Varivent N pipe DN65-162 PN25. Unit of measurement mm (in)

Tri-Clamp

NA Connect ISO2852 DN25-38 (1...1 1/2"), DIN32676 DN25-40


- Material: 316L
- Pressure rating: ≤ 25 bar (363 psi)
- Temperature: ≤ 150 °C (302 °F)
- Weight: 0.3 kg (0.66 lb)

NA Connect ISO2852 DN40-51 (2"), DIN32676 DN50

- Material: 316L
- Pressure rating: ≤ 25 bar (363 psi)
- Temperature: ≤ 150 °C (302 °F)
- Weight: 0.3 kg (0.66 lb)



A seal is not included in the delivery. The maximum temperature and maximum pressure depend on the clamping ring and sealing element used (according to the process connection design). The lowest value applies in each case.

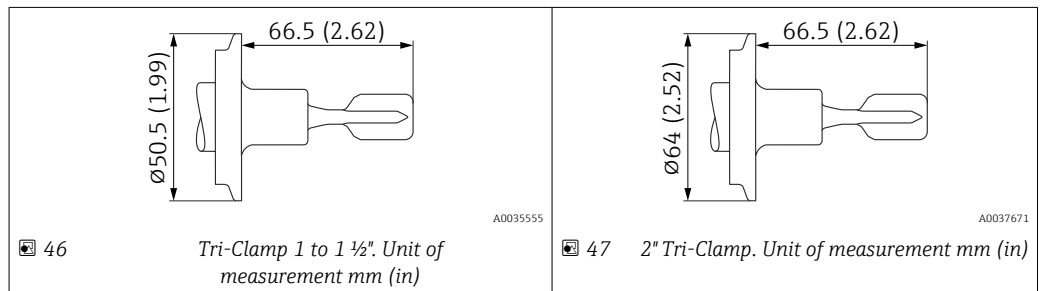

46

Tri-Clamp 1 to 1 1/2". Unit of measurement mm (in)

47

2" Tri-Clamp. Unit of measurement mm (in)

Weight

Basic weight: 0.65 kg (1.43 lb)

The basic weight comprises:

- Probe design: compact version
- Electronic insert
- Housing: single compartment, plastic with cover
- Thread, G 3/4



Differences in weight are caused by the housing and cover selected.

Housing

- Single compartment, aluminum, coated: 0.8 kg (1.76 lb)
- Single compartment; 316L, hygienic: 0.45 kg (0.99 lb)
- Dual compartment, L-shaped; aluminum, coated: 1.22 kg (2.69 lb)

Temperature spacer

0.6 kg (1.32 lb)

Pressure-tight feedthrough

0.7 kg (1.54 lb)

Pipe extension

- 1000 mm: 0.9 kg (1.98 lb)
- 50 in: 1.15 kg (2.54 lb)

Process connection

See "Process connections" section

Protective cover, plastic

0.2 kg (0.44 lb)

Protective cover, 316L

0.93 kg (2.05 lb)

Materials**Materials in contact with process***Process connection and pipe extension*

316L (1.4404 or 1.4435)

Tuning fork

316L (1.4435)

Seals

Scope of delivery including seal

- Ingot fitting, seal material: EPDM (in compliance with FDA, USP Class VI)
- Flush-mounted tank connection for installation in welding neck, seal material: silicone

Materials not in contact with process*Plastic housing*

- Housing: PBT/PC
- Dummy cover: PBT/PC
- Transparent cover: PA12
- Cover with sight glass: PBT/PC and PC
- Cover seal: EPDM
- Potential equalization: 316L
- Seal under potential equalization: EPDM
- Plug: PBT-GF30-FR
- M20 cable gland: PA
- Seal on plug and cable gland: EPDM
- Threaded adapter as substitute for cable glands: PA66-GF30
- Nameplate: plastic foil
- TAG plate: plastic foil, metal or provided by customer

Aluminum housing, coated

- Housing: aluminum EN AC 43400
- Dummy cover: aluminum EN AC 43400
- Cover seal materials: HNBR
- Cover seal materials: FVMQ
- Plug: aluminum
 - Plastic (PBT-GF30-FR) in Ex-free, Ex i or IS combination with cable gland, plastic, M20 thread or G ½ thread
- Nameplate: plastic foil
- TAG plate: plastic foil, stainless steel or provided by the customer
- M20 cable glands: select material (stainless steel, nickel-plated brass, polyamide)

Stainless steel housing, 316L, hygienic

- Housing: stainless steel AISI 316L (1.4404)
- Dummy cover: stainless steel AISI 316L (1.4404)
- Cover with polycarbonate sight glass optionally available. For dust ignition-proof applications, the sight glass is made of borosilicate.
- Cover seal materials: VMQ
- Plug: stainless steel or plastic
 - Plastic (PBT-GF30-FR) in Ex-free, Ex i or IS combination with cable gland, plastic, M20 thread or G ½ thread
 - Stainless steel for cable glands made of stainless steel or nickel or for Ex t, Ex ia IIIC
- Nameplate: stainless steel housing labeled directly
- TAG plate: plastic foil, stainless steel or provided by the customer
- M20 cable glands: select material (stainless steel, nickel-plated brass, polyamide)

Surface roughness	Surface roughness of the surface in contact with the process: Ra < 1.5 µm (59 µin) Optional: Ra < 0.3 µm (12 µin) mechanically polished (3-A, EHEDG)
-------------------	---

Mechanical construction of Density Computer FML621

Terminal	Pluggable screw terminals - power supply terminal is coded. The terminal range - 1.5 mm² (16 AWG) solid, 1 mm² (18 AWG) flexible with ferrules - applies to all connections.
----------	--

Dimensions	48 Housing for DIN rail as per IEC 60715. Unit of measurement mm (in)
------------	--


Slots with extension cards	49 Device with additional extension cards 1 Slot A, extension card (already included in the basic unit) 2 Slot B, extension card (optional or available as an accessory) 3 Slot C, extension card (optional or available as an accessory) 4 Slot D, extension card (optional or available as an accessory) 5 Slot E, extension card (already included in the basic unit)
----------------------------	---

Weight	Basic unit: 500 g (17.6 oz) Weight with all additional extension cards. Remote operating unit: 300 g (10.6 oz).
Materials	Housing: Polycarbonate plastic, UL 94V0

User interface of Density Computer FML621


-  An operating and display unit can also be used for commissioning the Density Computer FML621
- The operating and display unit can also be used for multiple devices
- An operating and display unit is absolutely essential for field adjustment

Display elements	Display 160x80 dot-matrix LCD with blue background lighting. In the event of an error, the color of the background lighting changes to red. It is possible to configure the background color. LED status indicator - Operation: 1 x green 2 mm (0.08 in) - Fault message: 1 x red 2 mm (0.08 in) Operating and display unit - optional or as an accessory - In addition, a display and operating unit with the following dimensions can be connected to the device in the panel-mounted housing: - W: 144 mm (5.67 in) - H: 72 mm (2.83 in) - D: 43 mm (1.69 in) - Connection to the integrated RS484 interface is performed using the connection cable (length = 3 m (9.84 ft)), which is included in the accessory kit - It is possible to operate the operating and display unit in parallel with a device-internal display in the FML621
-------------------------	--

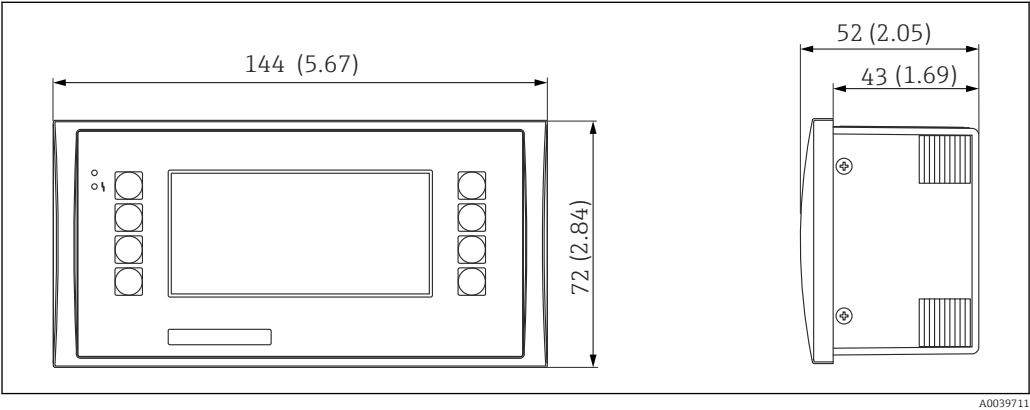

 50 Operating and display unit for panel mounting. Unit of measurement mm (in)

A0039711

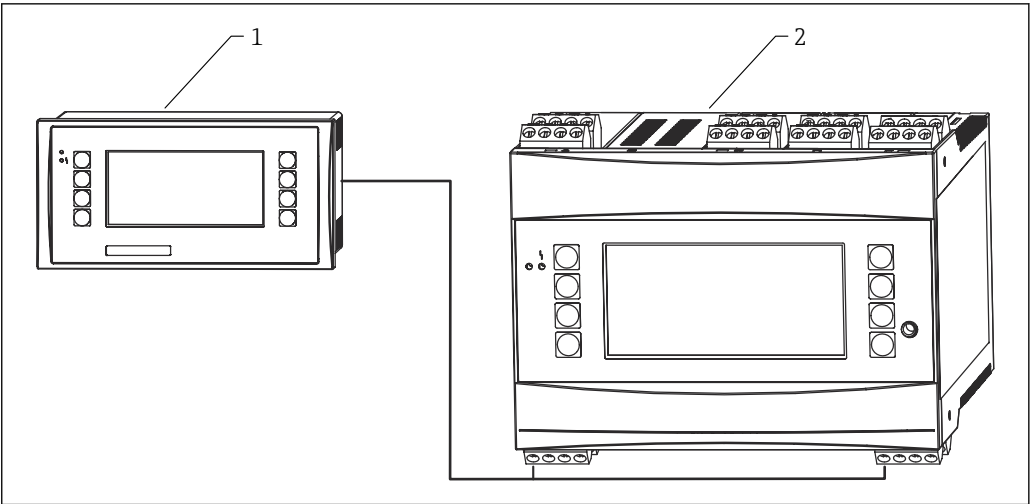

A0039717

51 Operating and display unit in panel-mounted housing

- 1 Operating and display unit
2 Basic unit

Operating elements	Eight soft keys at the front are used to interact with the display. The functions of the keys are shown on the display.
Remote operation	■ RS232 interface via mini jack plug socket 3.5 mm (0.14 in), configuration using a PC and ReadWin® 2000 PC software ■ RS485 interface
Real-time clock	■ Deviation: 30 min per year ■ Power reserve: 14 days

Certificates and approvals

CE mark	The measuring system meets the legal requirements of the applicable EU Directives. These are listed in the corresponding EU Declaration of Conformity along with the standards applied. Endress+Hauser confirms that the device has been successfully tested by applying the CE mark.
Ex approval	For available Ex approvals, see Product Configurator. All explosion protection data are listed in a separate document, which is available on request.
Other standards and guidelines	IEC 60529 Degrees of protection provided by enclosures (IP code) IEC 61010 Safety requirements for electrical equipment for measurement, control and laboratory use EN 61326 series EMC product family standard for electrical equipment for measurement, control and laboratory use NAMUR User association of automation technology in process industries
Additional information on Liquiphant Density	Material compliance for contact with food The device has been developed for food contact applications. Versions can be selected that meet the following requirements: ■ EU Food Contact Material (EC) 1935/2004 ■ US Food Contact Material FDA CFR 21 ■ CN Food Contact Material GB 4806

Hygienic design requirements

- Notes on installation and certification in accordance with 3-A and EHEDG:
 -  SD02503F document "Hygienic approvals"
- Information on 3-A and EHEDG-certified adapters:
 -  TI00426F document "Weld-in adapters, process adapters and flanges"
- 3-A and EHEDG-certified versions of the sensor are suitable for Cleaning-in-Place (CIP) and Sterilization-in-Place (SIP) without removing them from the plant. This means that the sensor does not need to be removed during cleaning. The maximum permitted pressure and temperature values for sensor and adapter must not be exceeded (see notes in this TI).
- ASME BPE

Compliance with requirements derived from cGMP

cGMP is suitable for wetted parts:

- The certificate is only available in English
- Materials of construction
- ADI-free based upon EMA/410/01 Rev.3 (TSE/BSE compliant)
- Polishing and surface finish
- Material/compound compliance table: USP, FDA

General material compliance

Endress+Hauser guarantees compliance with all relevant laws and regulations, including the current guidelines for materials and substances.

Examples:

- RoHS
- China RoHS
- REACH
- POP VO (Stockholm Convention)

For further information and general declarations of compliance, see the Endress+Hauser website www.endress.com

TSE (BSE) compliance (ADI free - Animal Derived Ingredients)

As the manufacturer, Endress+Hauser states:

- That the parts of this product in contact with the process are not made from materials derived from animals **or**
- at least comply with the requirements of guidelines outlined in EMA/410/01 rev. 3 (TSE (BSE) compliance).

Ordering information

Detailed ordering information is available from your nearest sales organization www.addresses.endress.com or in the Product Configurator at www.endress.com:

1. Select the product using the filters and search field.
2. Open the product page.
3. Select **Configuration**.



Product Configurator - the tool for individual product configuration

- Up-to-the-minute configuration data
- Depending on the device: Direct input of measuring point-specific information such as measuring range or operating language
- Automatic verification of exclusion criteria
- Automatic creation of the order code and its breakdown in PDF or Excel output format
- Ability to order directly in the Endress+Hauser Online Shop

TAG

Measuring point (TAG)

The device can be ordered with a tag name.

Location of tag name

Select in the additional specification:

- Stainless steel tag plate
- Paper adhesive label
- Tag provided by the customer
- RFID tag
- RFID tag + stainless steel tag plate
- RFID tag + paper adhesive label
- RFID tag + tag provided by the customer
- IEC 61406 stainless steel tag
- IEC 61406 stainless steel tag + NFC tag
- IEC 61406 stainless steel tag, stainless steel tag
- IEC 61406 stainless steel tag + NFC, stainless steel tag
- IEC 61406 stainless steel tag, plate supplied
- IEC 61406 stainless steel tag + NFC, plate supplied

Definition of tag name

Specify in the additional specification:

3 lines of maximum 18 characters each

The specified tag name appears on the selected plate and/or on the RFID tag.

Test reports, declarations and inspection certificates

All test reports, declarations and inspection certificates are provided electronically in the *Device Viewer*:

Enter the serial number from the nameplate (www.endress.com/deviceviewer)




Product documentation on paper

Test reports, declarations and inspection certificates in hard copy can optionally be ordered with feature 570 "Service", Version I7 "Product documentation on paper". The documents are then provided with the device upon delivery.

Accessories for Liquiphant Density

Protective cover for aluminum dual compartment housing


- Material: stainless steel 316L
- Order number: 71438303

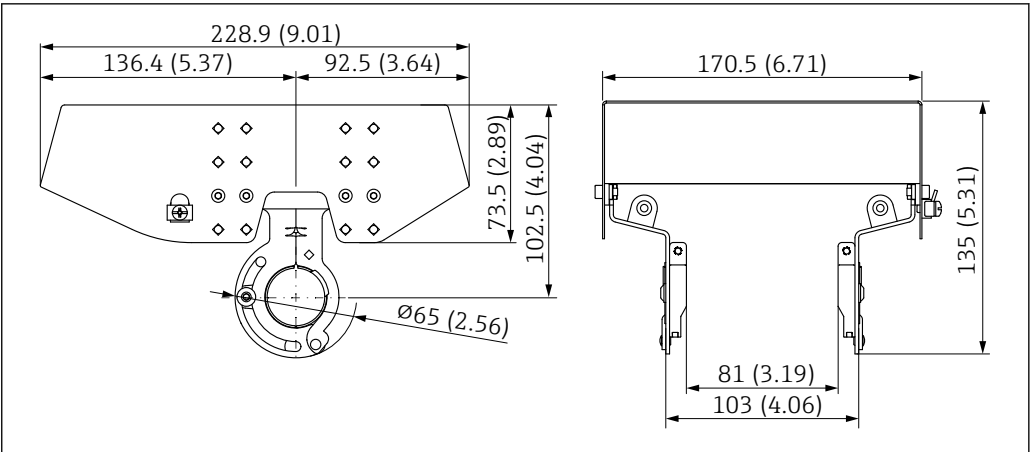

52 Protective cover for aluminum dual compartment housing. Unit of measurement mm (in)

Protective cover for aluminum single compartment housing

- Material: plastic
- Order number: 71438291

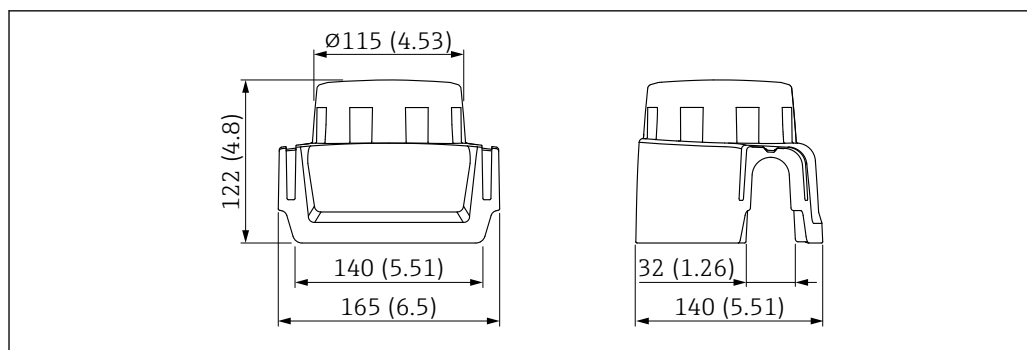


A0038280

53 Protective cover for aluminum single compartment housing. Unit of measurement mm (in)

Weld-in adapter

Various weld-in adapters are available for installation in vessels or pipes. The adapters are optionally available with inspection certificate 3.1 EN10204.

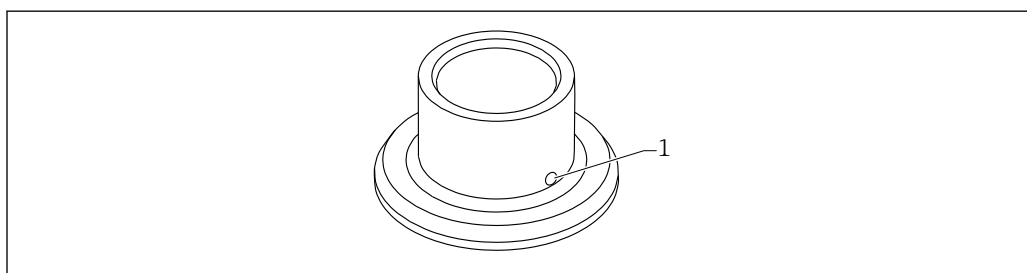

A0023557

54 Weld-in adapter with leakage hole (sample view)

1 Leakage hole

Weld in the weld-in adapter in such a way that the leakage hole is pointing downwards. This enables any leaks to be detected quickly.

- G 1, Ø53 mounting on the pipe
- G 1, Ø60 flush mount on the vessel
- G ¾, Ø55 flush mount
- G 1 sensor adjustable
- RD52 sensor adjustable



For detailed information, see "Technical Information" TI00426F (Weld-in adapters, process adapters and flanges)

Available in the Download Area of the Endress+Hauser website (www.endress.com/downloads).

M12 socket



The M12 sockets listed are suitable for use in the temperature range -25 to +70 °C (-13 to +158 °F).

M12 socket IP69

- Terminated at one end
- Angled
- 5 m (16 ft) PVC cable (orange)
- Slotted nut 316L (1.4435)
- Body: PVC
- Order number: 52024216

M12 socket IP67

- Angled
- 5 m (16 ft) PVC cable (gray)
- Slotted nut Cu Sn/Ni
- Body: PUR
- Order number: 52010285

Accessories for Density Computer FML621

General

RXU10-A1

Cable set for Density Computer FML621 for connecting to a PC or modem

FML621A-AA

Remote display for panel mounting:

- B: 144 mm (5.67 in)
- H: 72 mm (2.83 in)
- T: 43 mm (1.69 in)

RMS621A-P1

PROFIBUS interface

51004148

Adhesive label, printed, max. 2 x 16 characters

51002393

Metal sign for TAG number

51010487

Sign, paper, TAG 3 x 16 characters

Extension cards

The device can be extended with a maximum of three universal or digital or current or Pt100 cards.

FML621A-DA

Digital

- 6 x digital input
- 6 x relay output
- Kit with terminals and fixing frame

FML621A-DB

Digital, ATEX-approved

- 6 x digital input
- 6 x relay output
- Kit with terminals

FML621A-CA

2x U, I, TC

- 2x 0 to 20 mA or 4 to 20 mA per pulse
- 2x digital
- 2x SPST relay

FML621A-CB

Multifunction, 2x U, I, TC ATEX

- 2x 0 to 20 mA or 4 to 20 mA per pulse
- 2x digital
- 2x SPST relay

FML621A-TA

Temperature (Pt100/Pt500/Pt1000)

Complete, including terminals and securing frame

FML621A-TB

Temperature, ATEX-approved (Pt100/PT500/PT1000)

Complete, including terminals

FML621A-UA

Universal - PFM or pulse or analog or transmitter power supply

Complete, including terminals and securing frame

FML621A-UB

Universal ATEX-approved - PFM or pulse or analog or transmitter power supply

Complete, including terminals

PROFINET® Interface

Order code RMS621A-P2

Documentation



For an overview of the scope of the associated Technical Documentation, refer to the following:

- *Device Viewer* (www.endress.com/deviceviewer): Enter the serial number from the nameplate
- *Endress+Hauser Operations app*: Enter serial number from nameplate or scan matrix code on nameplate.

Standard documentation

Document type: Operating Instructions (BA)

Installation and initial commissioning – contains all the functions in the operating menu that are needed for a normal measuring task. Functions beyond this scope are not included.

Document type: Brief Operating Instructions (KA)

Quick guide to the first measured value – includes all essential information from incoming acceptance to electrical connection.

Document type: Safety Instructions, certificates

Depending on the approval, Safety Instructions are also supplied with the device, e.g. XA. This documentation is an integral part of the Operating Instructions.

The nameplate indicates the Safety Instructions (XA) that are relevant to the device.

Supplementary device-dependent documentation

Operating Instructions

BA00335F: Density Computer FML621

Special Documentation

- SD01622P: Weld-in adapter (installation instructions)
- TI00426F: Weld-in adapters, process adapters and flanges (overview)



71627519

www.addresses.endress.com

For more Endress+Hauser product please visit: www.ainstru.com

Mail: Sales@ainstru.com